



坚固的设计满足高要求的最终细度

锤式粉碎机适用于中硬、硬、脆、韧和纤维样品的破碎、粉碎和分散。

???RETSCH?????HM 200?????100 mm?????????????????
????????????0.8 mm??? HM200????????????????????????????
??HM 200????????????
???




[点击观看视频](#)

锤式破碎机 HM 200

高通量 & 明确的最终细度

?????????HM 200?????????HM 200?????????HM 200?????????125????????? SR 300?
SK 300?????????25 mm??HM 200??HM 200?????????100 mm?????????0.05 mm?0.1 mm?????????

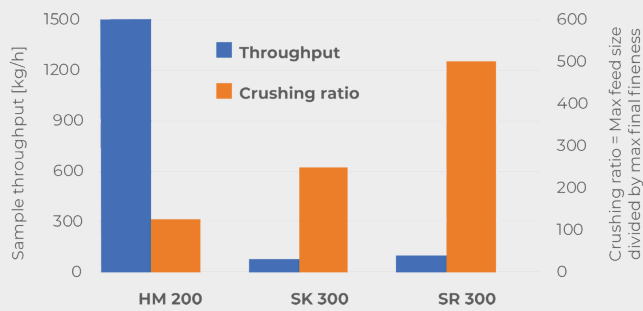

- | 优异的破碎性能
- | 高达1500公斤/小时的超高处理量
- | 固定转速3000 min⁻¹
- | 底筛孔径范围2 - 40 mm
- | 进样尺寸高达100 mm, 研磨细度 < 0.8 mm
- | 易于进入破碎室, 便于清洁
- | 适用于批量或连续研磨
- | 吸尘接头

锤式破碎机 HM 200

作用原理

HM 200 装有一个水平旋转轴，轴上安装有 10 个可自由摆动的锤子。这些锤子在研磨室内高速粉碎物料。与固定锤头相比，摆动锤头可最大限度地降低堵塞机器和磨损的风险。材料在被每个锤子击打和在研磨腔内弹跳时受到冲击而粉碎。一旦样品细到足以通过筛网，就会通过重力出口落入收集容器。



[点击观看视频](#)

锤式破碎机 HM 200

技术参数

应用	coarse and pre-crushing, desagglomeration
应用领域	food / feed, 农业, 化学 / 合成材料, 地质 / 冶金, 工程/电子, 建筑原料, 煤炭, 环境 / 资源回收利用, 玻璃/ 陶瓷
样品特征	中硬性, 硬的, 脆性的, 硬的, 含纤维的
粉碎原理:	碰撞, 剪切
最大进样尺寸	< 100 mm
最终出料粒度 *	< 0.8 mm
批次加料量*	up to 30 l depending on collecting receptacle
网频 50赫兹 (60赫兹) 下的转速	3000 min ⁻¹ (3500 min ⁻¹)
处理量*:	1500 kg/h
旋翼圆周线速度	31.4 m/s
转刀直径	200 mm
转刀类型	10 swing hammer
干磨	是
湿磨	无
低温研磨	无
研磨套件材料:	hardened steel 1.0503 (hammers); steel 1.0330 (sieves)
筛网尺寸:	2 / 3 / 4 / 5 / 6 / 8 / 10 / 15 / 40 mm
加配吸尘器	是
安装装备型供货	是
接收槽容积	10 l, 30 l, continuous outlet
驱动	三相电机
驱动功率	1.5 kW
电源数据:	不同电压
电源接头:	三相
防护类型	IP 54
机体尺寸 (宽x高x纵深)	700 x 1570 x 750 mm
净重	~ 150 kg
标准	CE

受样品材料性质和仪器配置/设定的影响

www.retsch.cn/hm200

ORDER DATA

HAMMER MILL HM 200

**Hammer Mill incl. rotor with swing hammers, wearing plates and 10 l collecting receptacle
(please order hopper and bottom sieve separately)**

21.753.0001	HM 200	3~ 400 V, 50 Hz
21.753.0002	HM 200	3~ 200-220 V, 60 Hz

其余可选的电源配置机型价格相同

HOPPER

23.785.0031	Long-stock hopper
23.785.0032	Standard hopper

BOTTOM SIEVES HM 200, STEEL

03.647.0563	梯形筛孔	0.50毫米
03.647.0564	梯形筛孔	1.00毫米
03.647.0592	梯形筛孔	1.50毫米
03.647.0565	圆形筛孔	2.00毫米
03.647.0566	圆形筛孔	3.00毫米
03.647.0567	圆形筛孔	4.00毫米
03.647.0568	圆形筛孔	5.00毫米
03.647.0569	圆形筛孔	6.00毫米
03.647.0570	圆形筛孔	8.00毫米
03.647.0571	圆形筛孔	10.00毫米
03.647.0572	方形筛孔	15.00 mm
03.647.0573	方形筛孔	40.00毫米

CYCLONE FOR HM 200

**CYCLONE SET WITH HOLDER, CONNECTING ELEMENT, LID WITH CONNECTOR FOR VACCUUM CLEANER
AND SAMPLE OUTLET
(PLEASE ORDER CASSETTE WITH OUTLET SEPERATELY)**

22.935.0039	Cyclone set with collecting receptacle 30 l
03.183.0138	Cassette with outlet for use with cyclone
02.605.0420	Adjustable steel connection tubing (316L) for cyclone set
05.707.0208	FDA certified hose, 1.5 m
22.748.0017	Industrial vacuum cleaner Attix, 230 V, 50/60 Hz

ADDITIONAL ITEMS HM 200

23.003.0039	Collecting receptacle, 10 litres, stainless steel
23.003.0038	Collecting receptacle, 30 litres, stainless steel
23.003.0037	Outlet for continuous operation
22.786.0036	Conversion kit for continuous operation, incl. standard hopper, outlet for continuous operation, collecting receptacle, 30 liters, stainless steel
22.748.0017	Industrial vacuum cleaner, 230 V, 50/60 Hz, incl. adapter (other electrical versions and spare filter bags available upon request)
23.711.0035	Wear plates, hardened steel
22.002.0028	Rotor incl. swing hammers