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

Manual

Planetary Ball Mill PM400



Translation



Copyright

© Copyright by
Retsch GmbH
Retsch-Allee 1-5
42781 Haan
Germany

Table of Contents

1	Notes on the Manual	6
1.1	Explanations of the Safety Instructions	7
1.2	General Safety Instructions	8
1.3	Repairs	9
2	Confirmation	10
3	Packaging, Transport and Installation	11
3.1	Packaging	11
3.2	Transport	11
3.3	Temperature fluctuations and condensed water	11
3.4	Conditions for the Installation Site	12
3.5	Installation of the Device	12
3.6	Type Plate Description	12
3.7	Electrical Connection	14
3.8	Creating interface connection	14
3.9	Transport	15
3.10	Installation of the Device	17
3.11	Installation of the Device	18
4	Technical Data	19
4.1	Use of the Device for the Intended Purpose	19
4.2	Number of grinding stations	20
4.3	Receptacle Volume	20
4.4	Feed Grain Size	20
4.5	Rated Power	20
4.6	Electrical Connection	20
4.7	Emissions	21
4.7.1	Noise levels	21
4.7.2	Measurement conditions	21
4.8	Degree of Protection	21
4.9	Protective Equipment	21
4.10	Dimensions and Weight	22
4.11	Required Floor Space	22
5	Operating the Device	23
5.1	Views of the Instrument	23
5.2	Overview table of the parts of the device	24
5.3	Operating elements and displays	25
5.4	Overview Table of the Operating Elements and the Display	25
5.5	Opening the device	25
5.6	Closing the device	25
5.7	Emergency Unlocking	26
5.8	Inserting the Grinding Jar	27
5.8.1	Clamping bolt	28
5.8.2	Inserting the grinding jar	31
5.8.3	Inserting the clamping unit	32
5.8.4	Function of the locking sleeve	32
5.9	Releasing the grinding jar clamping mechanism	33
5.10	Opening the Clamping Mechanism with the Opening Aid for the Clamping Unit	33
5.11	Power failure during grinding	34
5.12	Mahlbehälterauswahl für unterschiedlichen Probenmaterialien	34
5.13	Sample quantity	35
5.13.1	Guides for material quantity and balls	35
5.14	Ultrafine grinding	35
5.15	Stacking the grinding jars	36
5.15.1	Stacking the 50 ml and 80 ml grinding jars	36
5.15.2	Stacking the 25 ml and 12 ml grinding jars	36

5.16	Handling grinding jars	37
5.16.1	Carrying and gripping	37
5.16.2	Anti-rotation device	37
5.16.3	Heating the grinding jars	37
5.17	Grinding jar identification	38
5.18	Grinding jar cleaning	38
5.18.1	Drying the grinding jars	38
5.19	Opening and closing of the grinding jars with safety locking device	39
5.20	Wet Grinding with Highly Flammable Materials	40
6	Display and operation	42
6.1	Symbols in the Display Unit	42
6.2	Display unit – operation of the device	43
6.2.1	Adjustment options using the display menu	43
6.3	Direct access to the language menu	44
6.4	Menu structure	45
6.5	Operating modes	45
6.5.1	Manual operation	45
6.5.2	Program 01 to 10	45
6.5.3	Basic settings	46
6.6	Manual Mode	46
6.6.1	Grinding time	46
6.6.2	Speed	46
6.6.3	Interval	46
6.6.4	Direction reversal	46
6.6.5	Pause time	46
6.6.6	Save parameters	47
6.6.7	Start in	47
6.7	Programme Mode	47
6.7.1	Change program	47
6.7.2	Delete program	48
6.8	Basic settings	48
6.8.1	Automatic opening	48
6.8.2	Language	48
6.8.3	Brightness	48
6.8.4	Date	48
6.8.5	Time	49
6.8.6	Acoustic warning signal	49
6.8.7	Service	49
6.8.7.1	Operating hours	49
6.8.7.2	Software version of display	49
6.8.7.3	Software version of the controller	49
6.8.7.4	Update software	49
6.8.7.4.1	Safety notice	49
7	Fault messages	50
8	Installing additional equipment	51
8.1	Adapter for glass vials	51
8.1.1	Speed limits	53
8.1.2	Energy input	53
9	Cleaning, Wear and Maintenance	54
9.1	Maintenance	54
9.1.1	Servicing the clamping unit	55
9.1.2	Wear to the bases of the clamping unit	55
9.1.3	Rubber washer on the pressure plate	56

9.1.4	Wear to latching bracket	56
10	Return for Service and Maintenance	57
11	Disposal.....	58
12	Index	59

1 Notes on the Manual

This operating manual is a technical guide on how to operate the device safely and it contains all the information required for the areas specified in the table of contents. This technical documentation is a reference and instruction manual. The individual chapters are complete in themselves.

Familiarity (of the respective target groups defined according to area) with the relevant chapters is a precondition for the safe and appropriate use of the device.

This operating manual does not contain any repair instructions. If faults arise or repairs are necessary, please contact your supplier or get in touch with Retsch GmbH directly.

Application technology information relating to samples to be processed is not included but can be read on the Internet on the respective device's page at www.retsch.com.

Changes

Subject to technical changes.

Copyright

Disclosure or reproduction of this documentation, use and disclosure of its contents are only permitted with the express permission of Retsch GmbH.

Infringements will result in damage compensation liability.

1.1 Explanations of the Safety Instructions

In this Operating Manual we give you the following safety warnings

Serious injury may result from failing to heed these safety warnings. We give you the following warnings and corresponding content.

 **WARNING**

Type of danger / personal injury

Source of danger

- Possible consequences if the dangers are not observed.
 - **Instructions on how the dangers are to be avoided.**
-

We also use the following signal word box in the text or in the instructions on action to be taken:

 **WARNING**

Moderate or mild injury may result from failing to heed these safety warnings. We give you the following warnings and corresponding content.

 **CAUTION**

Type of danger / personal injury

Source of danger

- Possible consequences if the dangers are not observed.
 - **Instructions on how the dangers are to be avoided.**
-

We also use the following signal word box in the text or in the instructions on action to be taken:

 **CAUTION**

In the event of possible **property damage** we inform you with the word “Instructions” and the corresponding content.

NOTICE

Nature of the property damage

Source of property damage

- Possible consequences if the instructions are not observed.
 - **Instructions on how the dangers are to be avoided.**
-

We also use the following signal word in the text or in the instructions on action to be taken:

NOTICE

1.2 General Safety Instructions

 **CAUTION****Read the Operating Manual**

Non-observance of these operating instructions

- The non-observance of these operating instructions can result in personal injuries.
- **Read the operating manual before using the device.**
- **We use the adjacent symbol to draw attention to the necessity of knowing the contents of this operating manual.**



Target group : All persons concerned with the machine in any form

This machine is a modern, high performance product from Retsch GmbH and complies with the state of the art. Operational safety is given if the machine is handled for the intended purpose and attention is given to this technical documentation.

You, as the owner/managing operator of the machine, must ensure that the people entrusted with working on the machine:

- have noted and understood all the regulations regarding safety,
- are familiar before starting work with all the operating instructions and specifications for the target group relevant for them,
- have easy access always to the technical documentation for this machine,
- and that new personnel before starting work on the machine are familiarised with the safe handling of the machine and its use for its intended purpose, either by verbal instructions from a competent person and/or by means of this technical documentation.

Improper operation can result in personal injuries and material damage. You are responsible for your own safety and that of your employees.

Make sure that no unauthorised person has access to the machine.

 **CAUTION****Changes to the machine**

- Changes to the machine may lead to personal injury.
- **Do not make any change to the machine and use spare parts and accessories that have been approved by Retsch exclusively.**

*NOTICE***Changes to the machine**

- The conformity declared by Retsch with the European Directives will lose its validity.
 - You lose all warranty claims.
 - **Do not make any change to the machine and use spare parts and accessories that have been approved by Retsch exclusively.**
-

1.3 Repairs

This operating manual does not contain any repair instructions. For your own safety, repairs may only be carried out by Retsch GmbH or an authorized representative or by Retsch service engineers.

In that case please inform:

The Retsch representative in your country
Your supplier
Retsch GmbH directly

Your Service Address:

2 Confirmation

This operating manual contains essential instructions for operating and maintaining the device which must be strictly observed. It is essential that they be read by the operator and by the qualified staff responsible for the device before the device is commissioned. This operating manual must be available and accessible at the place of use at all times.

The user of the device herewith confirms to the managing operator (owner) that (s)he has received sufficient instructions about the operation and maintenance of the system. The user has received the operating manual, has read and taken note of its contents and consequently has all the information required for safe operation and is sufficiently familiar with the device. As the owner/managing operator you should for your own protection have your employees confirm that they have received the instructions about the operation of the machine.

I have read and taken note of the contents of all chapters in this operating manual as well as all safety instructions and warnings.

User

Surname, first name (block letters)

Position in the company

Signature

Service technician or operator

Surname, first name (block letters)

Position in the company

Place, date and signature

3 Packaging, Transport and Installation

3.1 Packaging

The packaging has been adapted to the mode of transport. It complies with the generally applicable packaging guidelines.

NOTICE

Storage of packaging

- In the event of a complaint or return, your warranty claims may be endangered if the packaging is inadequate or the machine has not been secured correctly.
- **Please keep the packaging for the duration of the warranty period.**

3.2 Transport

NOTICE

Transport

- Mechanical or electronic components may be damaged.
- **The machine may not be knocked, shaken or thrown during transport.**

NOTICE

H0014

Complaints

- The forwarding agent and Retsch GmbH must be notified immediately in the event of transport damage. It is otherwise possible that subsequent complaints will not be recognised.
- **Notify your forwarding agent and Retsch GmbH within 24h**

3.3 Temperature fluctuations and condensed water

NOTICE

Temperature fluctuations

The machine may be subject to strong temperature fluctuations during transport (e.g. aircraft transport)

- The resultant condensed water may damage electronic components.
 - **Protect the machine from condensed water.**
-

3.4 Conditions for the Installation Site

Ambient temperature: 5°C to 40°C

NOTICE

Ambient temperature

- Electronic and mechanical components may be damaged and the performance data alter to an unknown extent.
- **Do not exceed or fall below the permitted temperature range of the machine (5°C to 40°C / ambient temperature).**

3.5 Installation of the Device

Installation height: maximum 2000 m above sea level

3.6 Type Plate Description

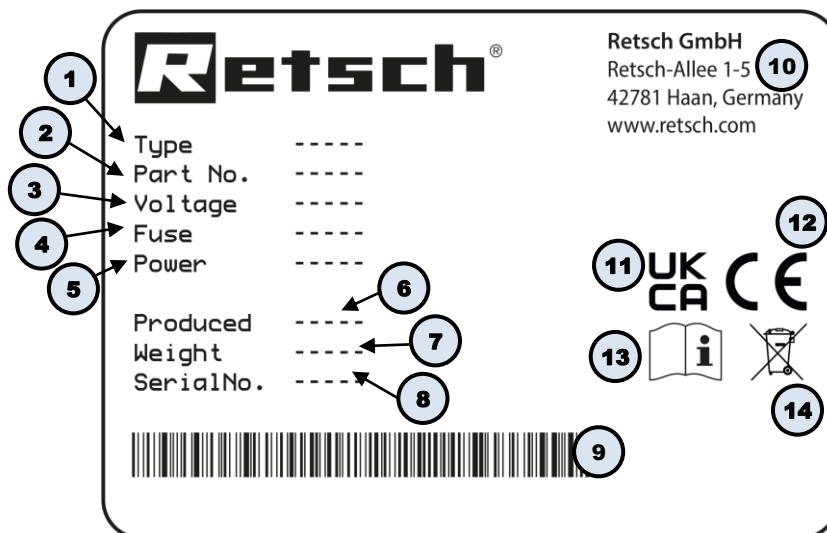


Fig. 1: Type plate

- 1 Device designation
- 2 Part number
- 3 Power version, Mains frequency
- 4 Fuse type and fuse strength
- 5 Capacity, Amperage
- 6 Year of production
- 7 Weight
- 8 Serial number
- 9 Bar code
- 10 Manufacturer's address
- 11 UKCA marking
- 12 CE marking
- 13 Safety warning: Read the manual
- 14 Disposal label

① In the case of queries please provide the device designation **(1)** or part number **(2)**, as well as the serial number **(8)** of the device.

3.7 Electrical Connection

⚠ WARNING

When connecting the power cable to the mains supply, use an external fuse that complies with the regulations applicable to the place of installation .

- Please check the type plate for details on the necessary voltage and frequency for the device.
- Make sure the levels agree with the existing mains power supply.
- Use the supplied connection cable to connect the device to the mains power supply.
- Make sure that the voltage and frequency of your mains connection corresponds to that on the type plate of the RS200.
 - The mains connection must be fused to at least 16A
 - An electrical connection without protective earth PE is not permitted.

The drive of the device is equipped with a frequency converter . In order to satisfy the EMC Directive , this is fitted with a mains filter and shielded cables to the motor. If your mains connection includes a residual current protection device, the suppressor capacitor wiring of the frequency converter when this is switched on (it is switched on by closing the grinding chamber hood) can lead to accidental triggering of the residual current protection device without any error being present on the device or in the mains installation.

In accordance with the state of the art, selective all current sensitive residual current protection devices are recommended for such cases. The tripping current must be sufficiently dimensioned because capacitive compensating current (shielded cable, mains filter) which only occurs for a short time can easily lead to accidental triggering.

In certain circumstances it may be necessary to operate the device without a residual current protection device. It is then necessary, however, to check that this does not contravene the local regulations of the electricity company or other institutions and the applicable standards.

3.8 Creating interface connection



Fig. 2: Serial interface

- When updating the device control or display software, a connection with the PC is established using the RS232 interface (**SC**).

NOTICE

The interface cables must not be longer than 2.5m. Longer cables can lead to faults in data transmission.

3.9 Transport

 **WARNING**

Serious personal injury

Falling loads

- The appliance is very heavy and can therefore cause serious personal injuries if it falls down.
- **Lifting above head height is not permissible!**

NOTICE

Transport

- Mechanical or electronic components may be damaged.
- **The machine may not be knocked, shaken or thrown during transport.**

NOTICE

N1.0018

Transportation lock

Transport without transportation lock, or operation with transportation lock

- Mechanical components may be damaged.
- **Only transport the device with mounted transportation lock.**
- **Do not operate the device with built-in transportation lock.**

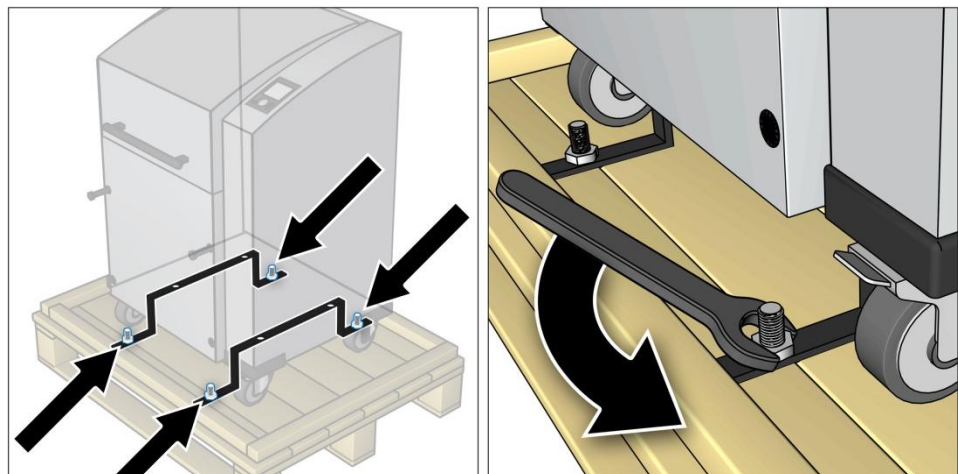


Fig. 3: Unscrewing the transport lock from the transport pallet

The device is secured to the transport pallet by the transport lock and four nuts.

- Use a 13mm spanner to unscrew the four nuts.

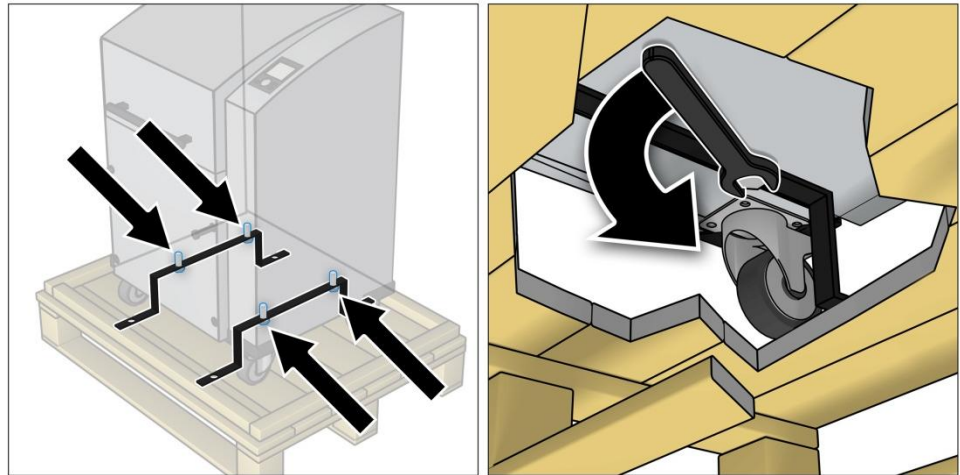


Fig. 4: Removing the transport lock from the device

Four bolts secure the transport lock underneath the device.

- Use a 13mm spanner to unscrew the four bolts.

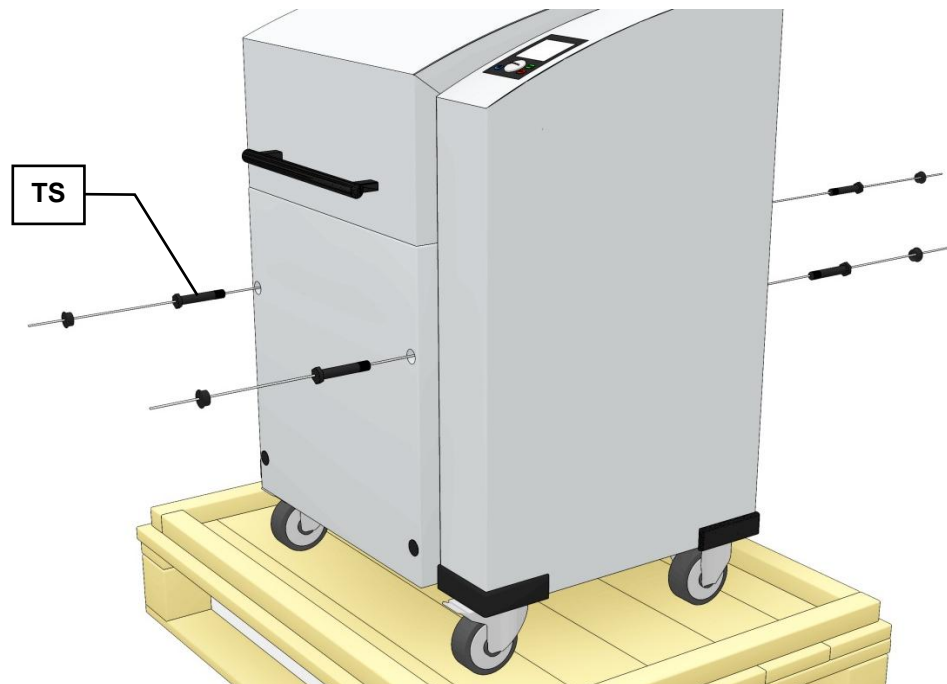


Fig. 5: Mounting the transport screws

The device should only be lifted and transported using the 4 transport screws (**TS**) provided.
Net weight approx. 210 kg

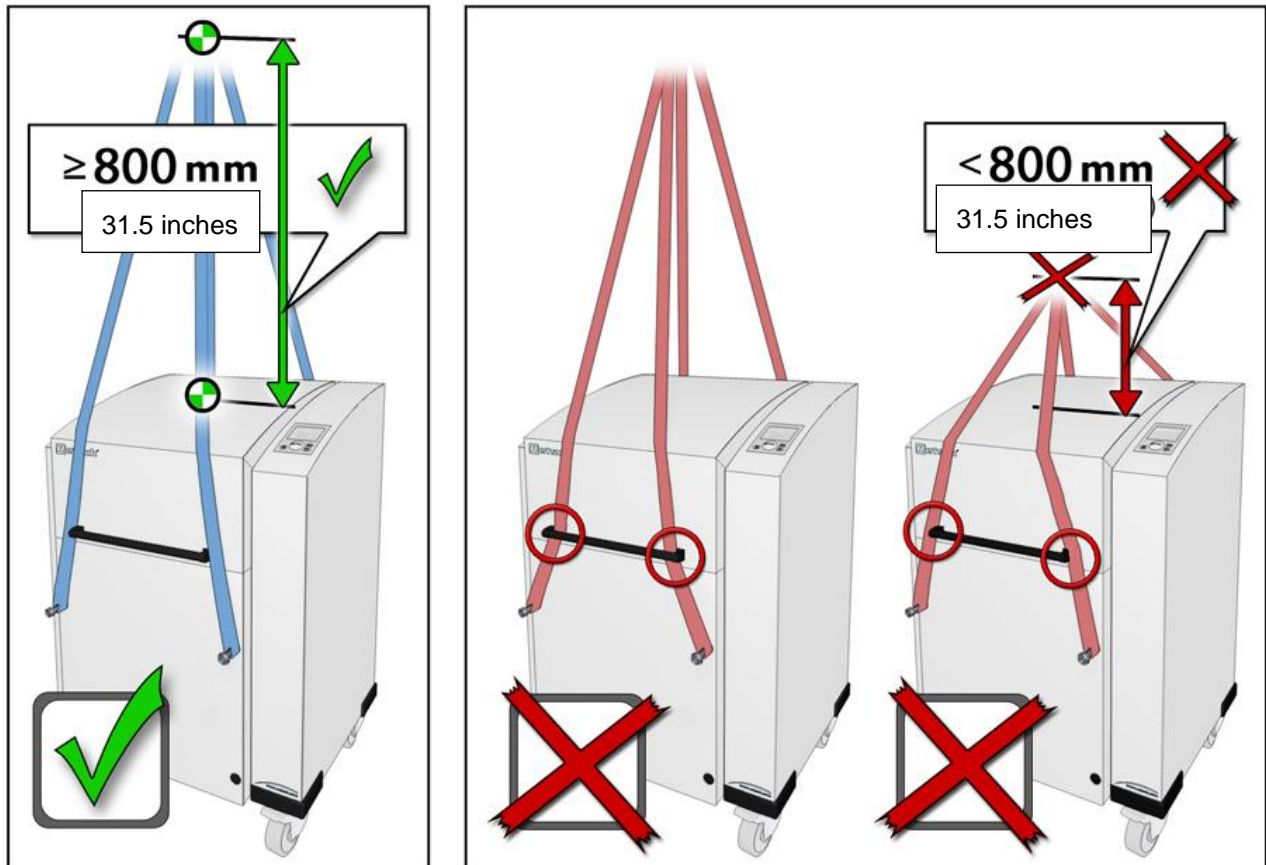


Fig. 6: Attaching the hoist

3.10 Installation of the Device

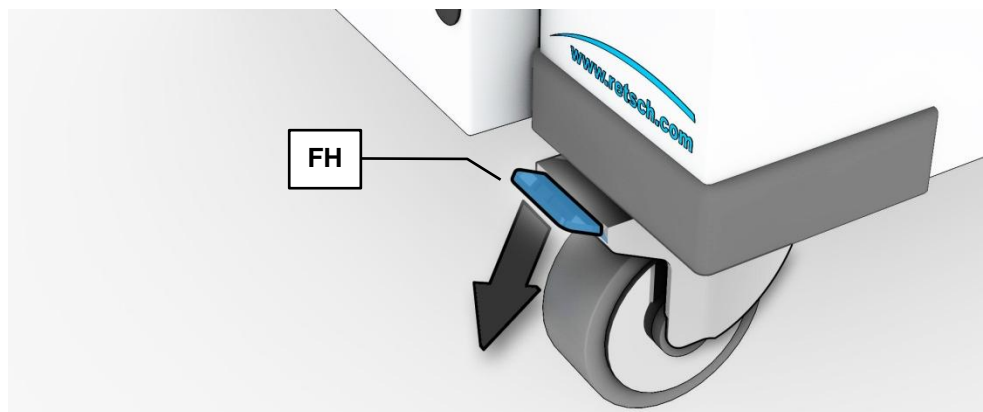


Fig. 7: Locking the transport rollers

- Place the device on a firm surface.
Please refer to the “Technical Data” chapter for further parameters.
The device must be locked before it is put into operation.
- Press the locking lever (FH) of the two front rollers down.

3.11 Installation of the Device

Installation height: maximum 2000 m above sea level

NOTE

Installation

- Depending on the operating status of the mill, there may be slight vibrations.
 - **Place the mill on an even, flat and balanced supporting surface only. The supporting surface must be stable and must not vibrate.**
-

NOTICE

Installation of the machine

- It must be possible to disconnect the machine from the mains at any time.
 - **Install the machine such that the connection for the mains cable is easily accessible.**
-

4 Technical Data

4.1 Use of the Device for the Intended Purpose

 **CAUTION**

1.V0004

Risk of explosion or fire

Changing sample properties

- Consider that the properties and therefore also the hazardousness of your sample can change during the grinding process.
- **Do not use any substances in this device which carry the risk of explosion or fire.**

 **CAUTION****Risk of explosion or fire**

– On account of its design, the device is not suitable for use in hazardous (potentially explosive) atmospheres.

- **Do not operate the device in a hazardous atmosphere.**

 **CAUTION****Danger of personal injury**

Dangerous nature of the sample

- Depending on the dangerous nature of your sample, take the necessary measures to rule out any danger to persons.
- **Observe the safety guidelines and datasheets of your sample material.**



Target group: Operating companies, operators

Machine type designation: PM400

Retsch ball mills are used to grind and mix soft, medium hard and extremely hard, brittle and fibrous materials. Dry and wet grinding are possible. Minerals, ores, alloys, chemicals, glass, ceramics, parts of plants, soil, sewage sludge, house or industrial waste and many other substances can be ground simply, quickly and without loss. The ball mills are used successfully in almost all areas of industry and research. This applies in particular where there are high demands in terms of hygiene, speed, fineness and reproducibility.

Only grinding jars in Comfort design (discontinued summer 2023) and EasyFit design are compatible with the device.

Grinding jars in the sizes 12 ml, 25 ml, 50 ml, 80 ml, 125 ml, 250 ml and 500 ml can be used.

Stacking of grinding jars is possible in the sizes 12 ml, 25 ml, 50 ml and 80 ml.

In order to insert and stack specific grinding jars, you may need the adapters which are available as accessories.

NOTICE

Area of use of the machine

- This machine is a laboratory machine designed for 8-hour single-shift operation.
 - **This machine may not be used as a production machine nor is it intended for continuous operation.**
-

4.2 Number of grinding stations

2 or 4 depending on the model

The grinding stations must be operated with the identical grinding jars and with the same weight during each grinding process.

NOTICE

H0068

Strong vibration and loud noise

Uneven load

- The device can produce very high levels of vibrations and noise when loading is uneven.
 - **Always use 2 or 4 grinding jars opposite each other.**
 - **When using 2 grinding jars, the two free grinding stations must be secured using a clamping bolt and a lead frame.**
 - **The grinding stations must be operated with the identical grinding jars and with the same weight during each grinding process.**
 - **Switch the device off immediately if there are high levels of vibration and noise, and check the number and net weight of the jars.**
-

4.3 Receptacle Volume

Up to 4 times 300ml, depending on the grinding jar volume.

4.4 Feed Grain Size

Up to <10 mm, however this depends on the material.

4.5 Rated Power

1500W

Make sure that the voltage and frequency of your mains connection correspond to that on the type plate of the device. The mains connection must be fused to at least 16A .

4.6 Electrical Connection

WARNING

When connecting the power cable to the mains supply, use an external fuse that complies with the regulations applicable to the place of installation .

- Please check the type plate for details on the necessary voltage and frequency for the device.

- Make sure the levels agree with the existing mains power supply.
- Use the supplied connection cable to connect the device to the mains power supply.

The drive for the appliance is equipped with a frequency converter. To fulfil the EMC directive, this is equipped with a mains filter and shielded cables to the drive. If your mains connection for the device includes a residual current device, the fault protection circuit on the frequency inverter may cause the residual current device to trip incorrectly when it is switched on without there being a fault in the device or in the mains installation.

According to the current state of the art, selective, all-current sensitive residual current devices of type B or B+ are recommended in such cases. The tripping current must be sufficiently dimensioned, as capacitive equalising currents that occur only briefly can easily lead to false tripping when switching the unit on.

4.7 Emissions



Damage to hearing

The level of noise can be high depending on the type of material, the knife used, the speed set and the duration of the grinding process.

- Noise that is excessive in terms of level and duration can cause impaired or permanently damaged hearing.
- **Ensure suitable sound-proofing measures or wear hearing protection.**



4.7.1 Noise levels

Noise measurement in accordance with DIN 45635-31-01-KL3

The noise levels are largely influenced by the machine speed, the grinding jar size and the diameter of the grinding balls used.

Workplace-related emissions L_{pAeq} = up to 85dB(A)

4.7.2 Measurement conditions

Grinding set: 4x 500ml special steel with 5 balls each Ø30mm TC (tungsten carbide)

Sample material: quartz, 135g each

Speed: 380 min⁻¹

4.8 Degree of Protection

IP40

4.9 Protective Equipment

This device is equipped with automatic lid closing which prevents it being started in an unsafe state.

- The device can only be started with closed lid.
- The lid can only be opened when the device has come to a halt.

4.10 Dimensions and Weight

Height: up to approx. 1220 mm / Width: 836 mm / Depth: up to approx. 780 mm
Weight : PM400 net approx. 290 kg

4.11 Required Floor Space

Height (open cover): 1900mm / Width: 1400mm / Depth: 900mm;

NOTICE

A clearance distance of 100mm is necessary at the back to allow the fans to operate.

5 Operating the Device

5.1 Views of the Instrument

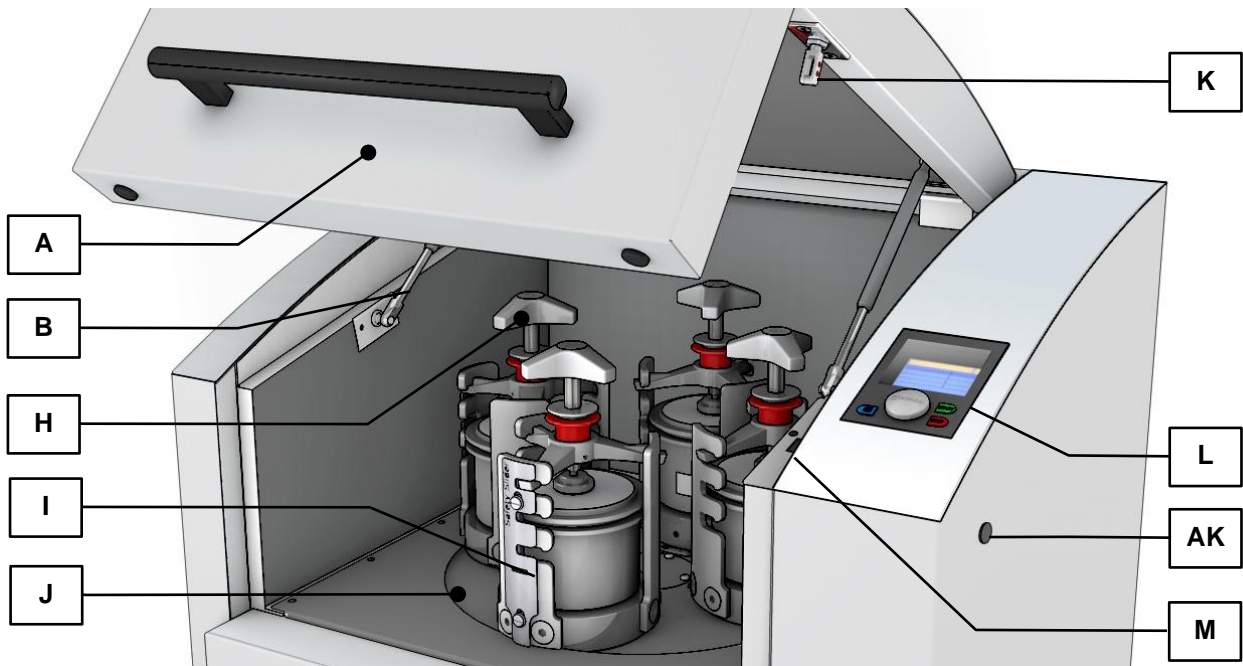


Fig. 8: View of the front of the device

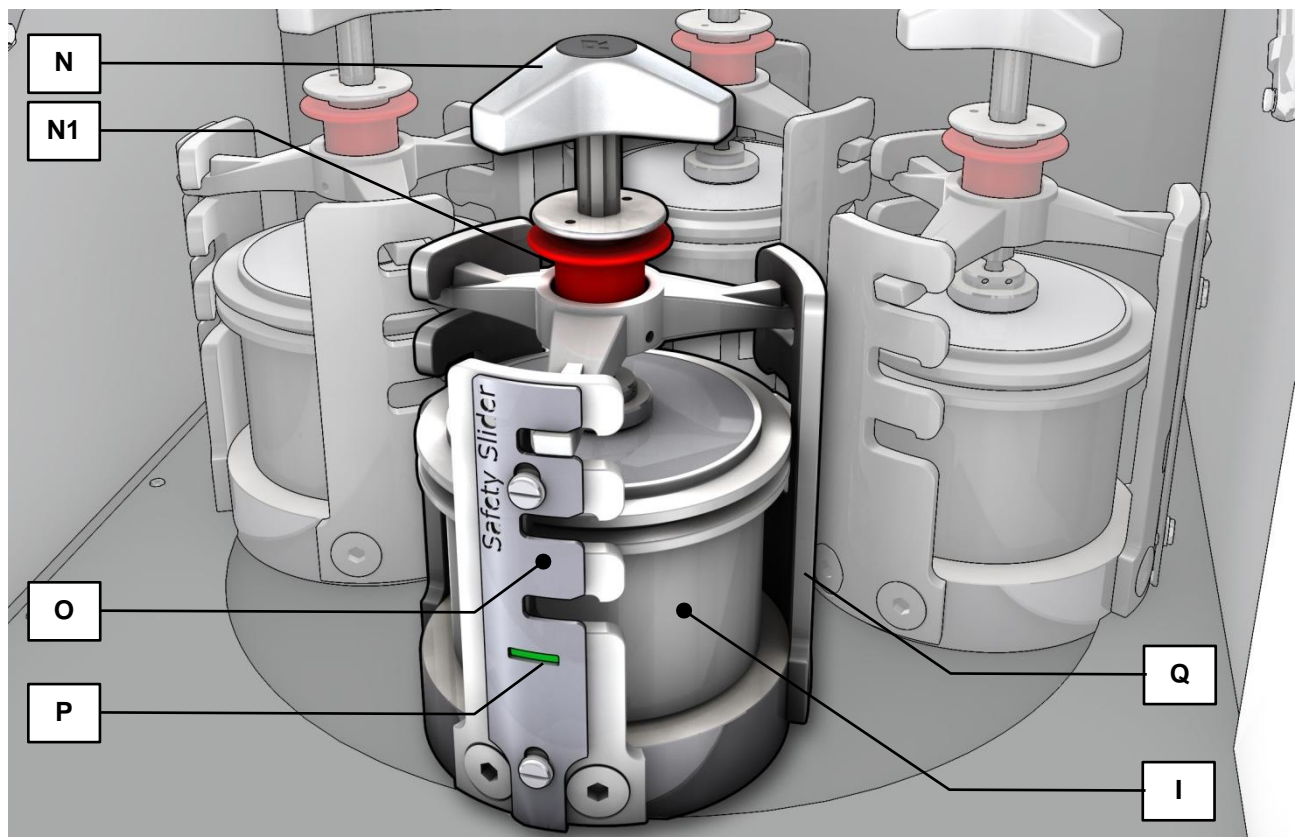


Fig. 9: Close-up of the grinding chamber

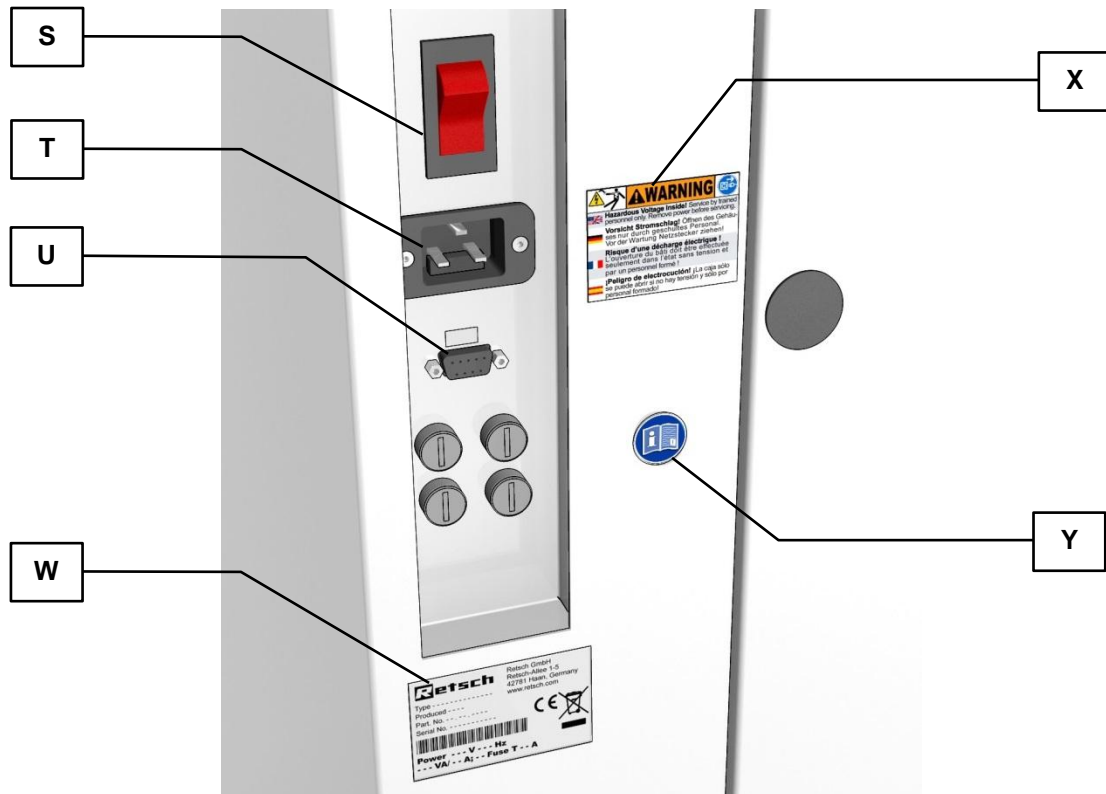


Fig. 10: View of the back of the device

5.2 Overview table of the parts of the device

Element	Description	Function
A	Cover	Closes the grinding chamber
B	Damper for cover	Secures the cover when it is open
H	Clamping unit for grinding chamber	Clamps grinding jar on the support
I	Grinding jar	Sample vessel
J	Turntable	Support for the grinding jar plate
K	Pin closing mechanism	Safety lock for the cover
L	Control panel	Device controller
M	Closing mechanism	Support for pin closing mechanism
N	Handle for clamping unit (lead frame)	Handle to clamp the grinding jars
N1	Locking sleeve	Prevents the twist grip from being turned accidentally
O	Safety Slider	Safety component – checking whether the clamping unit is inserted and clamped
P	Display window	Displays the safety function
Q	Latching bracket	Supports the clamping unit
S	Main switch	Disconnects the device from the mains
T	IEC socket	Power supply for the power cable
U	Serial interface	Interface for communication with the device
W	Type plate	Description of device parameters
X	Warning sign to pull out the plug	Warning of electric shock
Y	Sign with instruction to read the operating manual	Sign pointing out the need to read the operating manual

5.3 Operating elements and displays



Fig. 11: View of the control panel

5.4 Overview Table of the Operating Elements and the Display

Element	Description	Function
F	Control knob (rotary/push button)	Rotary control to operate the device settings
G	Button to open the cover	Unlocks the cover
C	Display	Displays the control functions and parameters
D	START button	Starts grinding
E	STOP button	Stops grinding

5.5 Opening the device

The following steps are necessary in order to be able to use and clamp the grinding jar.

- Connect the device to the mains.
- Switch on the main switch at the back of the device.

- Press the  button.

The safety lock opens and the lid can be lifted up

5.6 Closing the device

It is only possible to lock the grinding chamber if the device has been connected to the power supply and the main switch at the back of the device has been switched on.

- Shut the housing lid.
 - A sensor detects the closing pin of the housing lid and the motorised lid closing mechanism is switched on.
 - The housing lid is locked automatically.

5.7 Emergency Unlocking

CAUTION

Emergency Unlocking

Drive continuing to run

- There is a substantial risk of injury if the drive and associated device parts run on a long time without being braked!
- **Activate the emergency unlocking only when the machine has come to a complete stop and is disconnected from the power supply.**

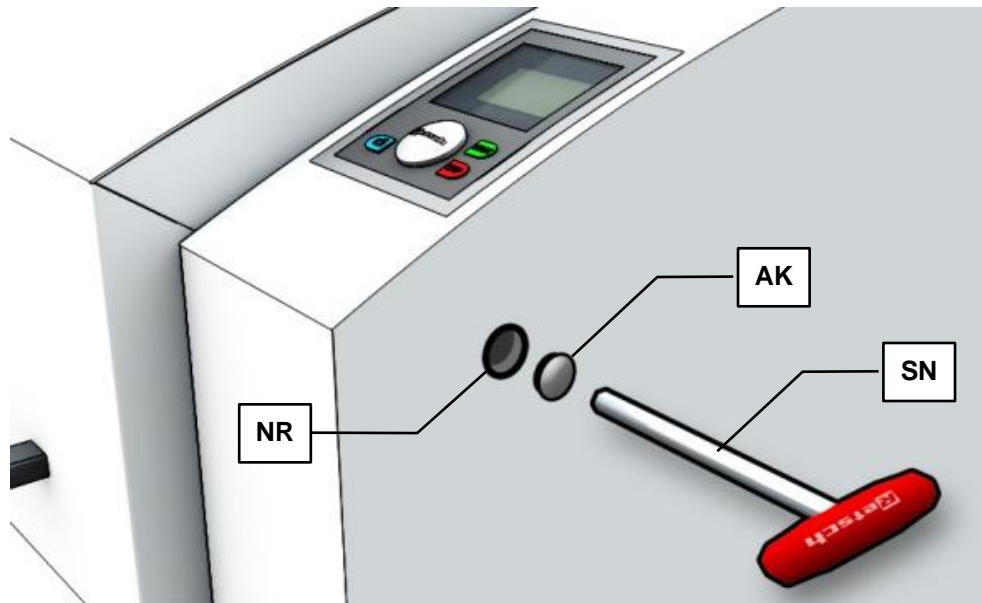


Fig. 12: Emergency release

A key is provided with device delivery. This can be used to open the device manually in the case of a power failure .

- Remove the cap (AK).
- Place the key (SN) in the opening (NR).
- To unlock the gear mechanism, it is necessary to push the key in further using a little force.

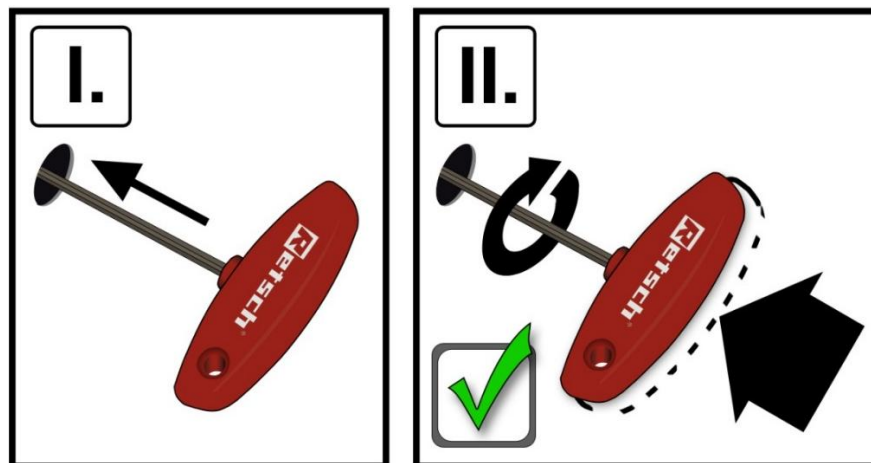


Fig. 13: Emergency release procedure

- At the same time as pushing the key (**SN**) in, rotate it in a clockwise direction as far as it will go.
- The lock is open and the lid can be lifted up.

5.8 Inserting the Grinding Jar

CAUTION

V0049

Ejected objects

Grinding jar not clamped

- Grinding jars or clamping mechanisms can be ejected. This results in a risk of injury.
- **Never leave loose clamping devices without clamped grinding jars in the grinding jar support.**
- **Before starting the machine, ensure that all grinding jars are clamped.**
- **Ensure that the red sleeve of the clamping mechanism is clicked into place.**

- **In the case of extended grinding processes, check that the grinding jars are secure at the following intervals:
After 3 min, after 1h, after 5h then every 10-12h.**



CAUTION

Scalding/burns

Hot grinding jar

- Depending on the grinding process, the material being ground and accordingly the grinding jar can become very hot.
- **Wear appropriate protection always when touching the grinding jar if it is hot.**
- **Never open hot grinding jars! Allow the grinding jars to cool down to room temperature before opening.**

NOTICE

H0068

Strong vibration and loud noise

Uneven load

- The device can produce very high levels of vibrations and noise when loading is uneven.
- **Always use 2 or 4 grinding jars opposite each other.**
- **When using 2 grinding jars, the two free grinding stations must be secured using a clamping bolt and a lead frame.**
- **The grinding stations must be operated with the identical grinding jars and with the same weight during each grinding process.**
- **Switch the device off immediately if there are high levels of vibration and noise, and check the number and net weight of the jars.**

NOTICE

H0011

Wear or damage to the grinding set

Different materials

- During the operation of the grinding set consisting of different materials, increased wear or damage to the grinding set is possible.
- **Only use grinding sets, where all parts are made of the same material.**

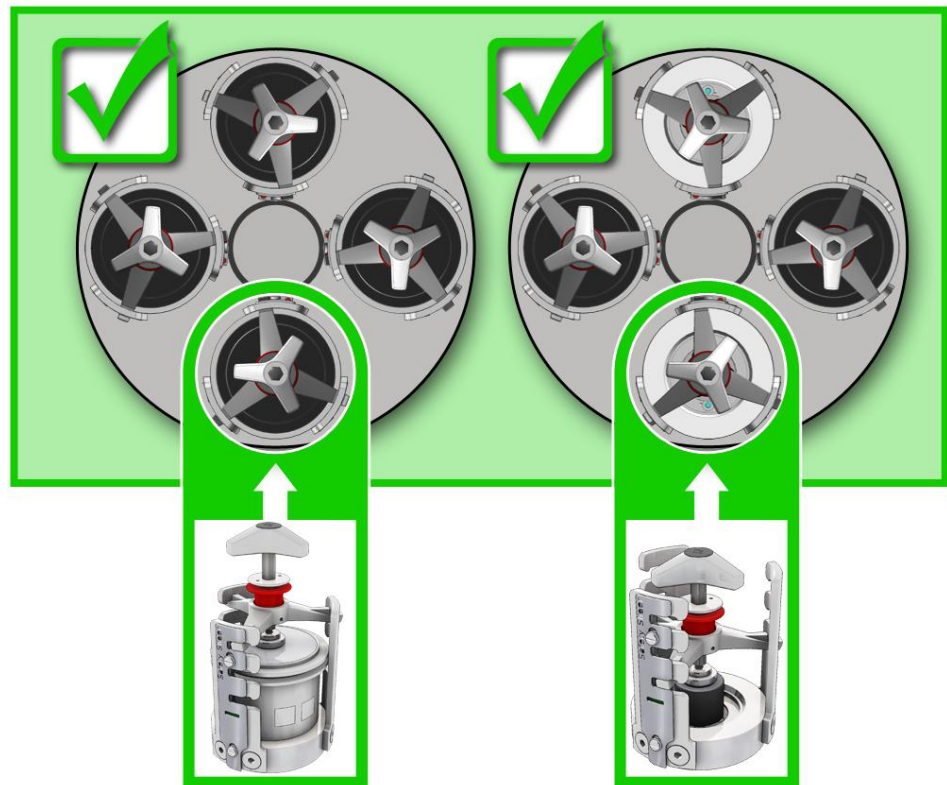
5.8.1 Clamping bolt

Fig. 14: Correct loading of the device – 4 or 2 grinding jars

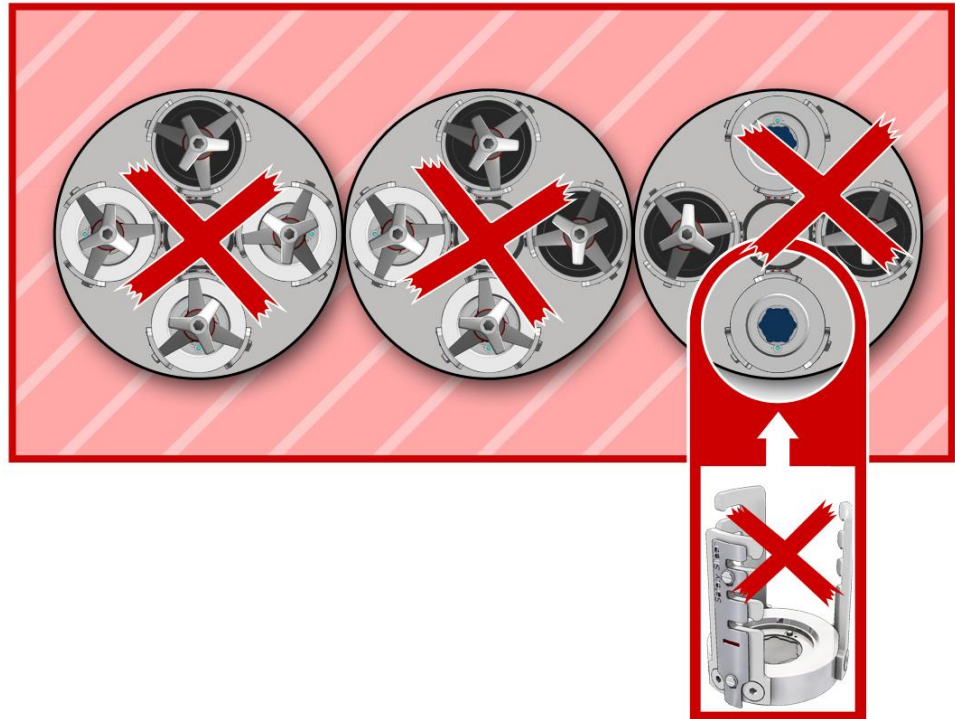


Fig. 15: Incorrect loading of the device

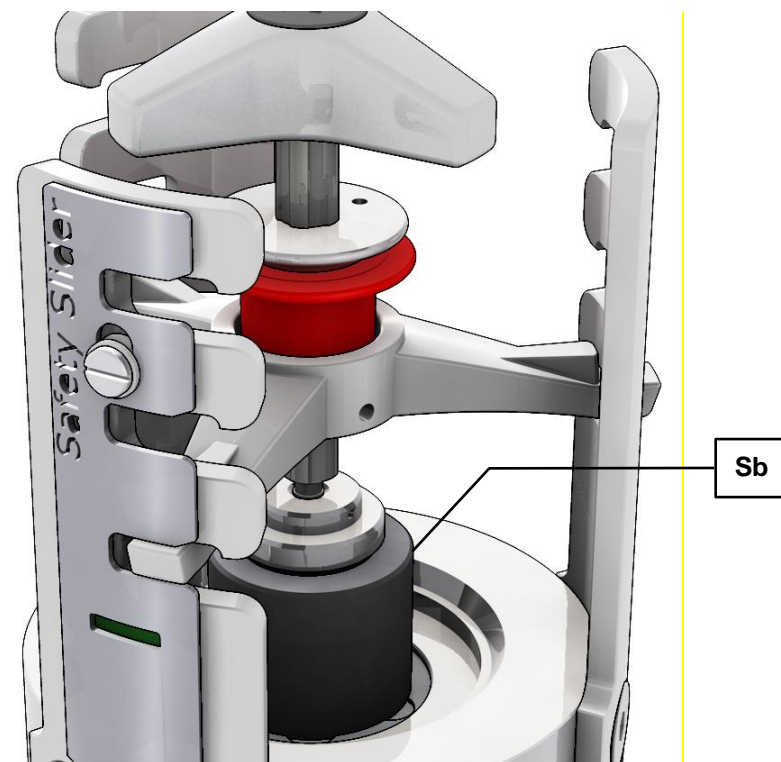


Fig. 16: Clamping bolt

When grinding with 2 grinding jars these must be inserted so they are opposite each other. The two remaining empty grinding stations must each be secured by a clamping bolt (**Sb**) and a clamping unit.

Otherwise the safety slider function will stop the device from starting.

Strong oscillation of the machine indicates poor weight balance or incorrectly clamped grinding jars!

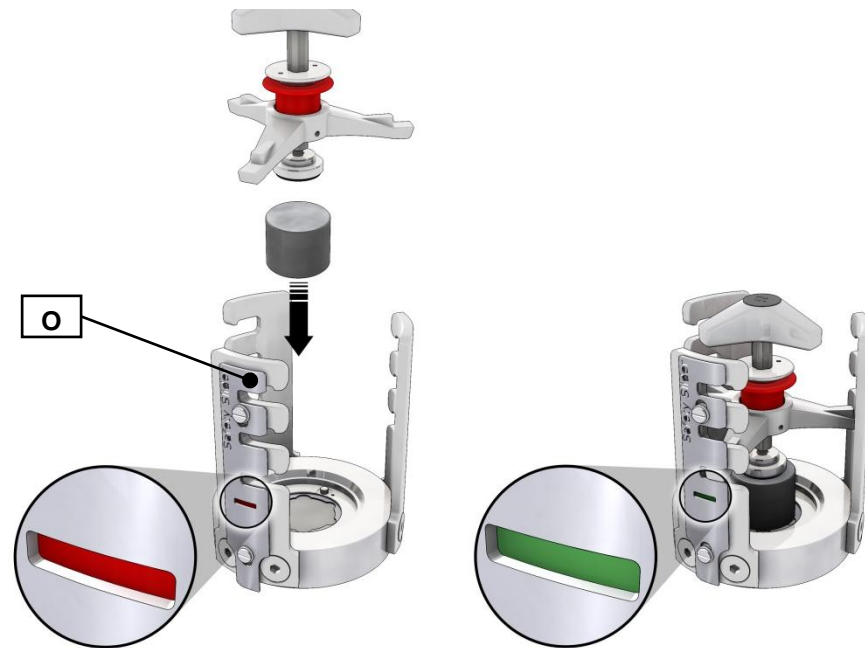


Fig. 17: Clamping bolt activates the Safety Slider

- When using two grinding jars, place a clamping bolt and a clamping unit in the two empty opposite grinding stations.
 - The insertion and clamping of the clamping unit pushes the Safety Slider (O) upwards, thereby activating it.
- After the machine has started, the device checks this safety function for approx. 15 seconds. The device aborts the starting process and emits an error message if
- you have not clamped the clamping unit;
 - you have not inserted any clamping unit or clamping bolts; or
 - if the grinding jar is missing.

5.8.2 Inserting the grinding jar

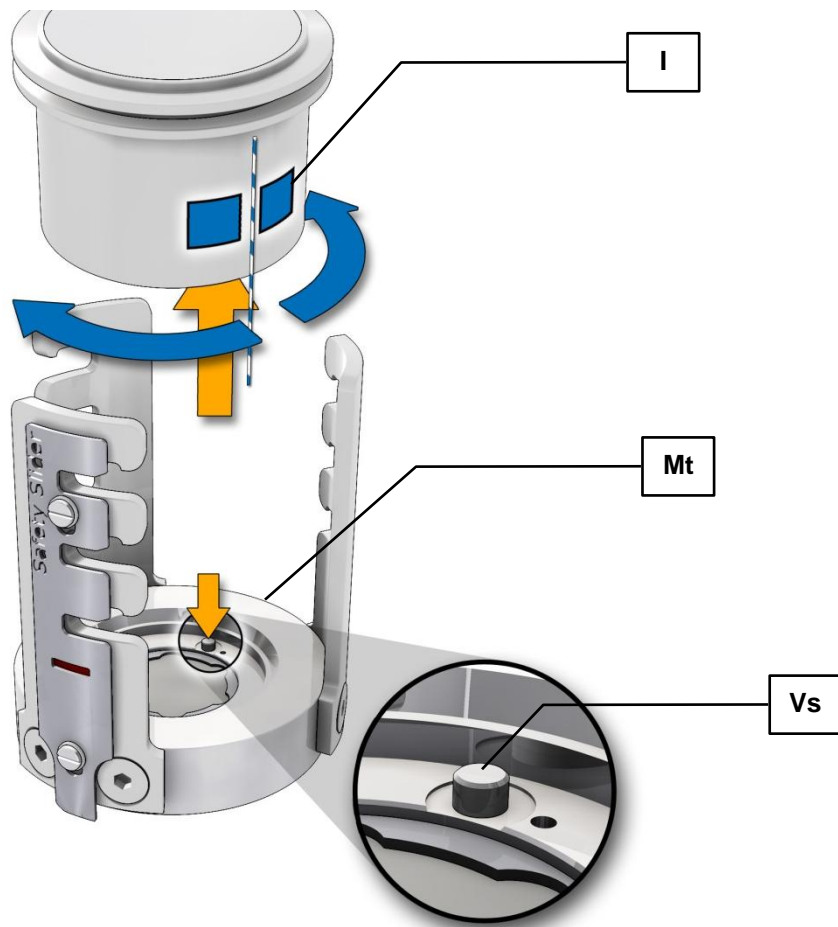


Fig. 18: Clamping the grinding jar

- If necessary clean the grinding jar base (**Mt**) and the anti-rotation pin (**Vs**).
- Turn the grinding jar label (**I**) to the same side as the anti-rotation pin.
- Place the grinding jar in the grinding jar support.
 - Pay attention to the anti-rotation device when using 250 and 500 ml grinding jars. The hole drilled in the bottom of the grinding jar to support the anti-rotation pin (**Vs**) is situated on the side of the lettering areas (**I**).

5.8.3 Inserting the clamping unit

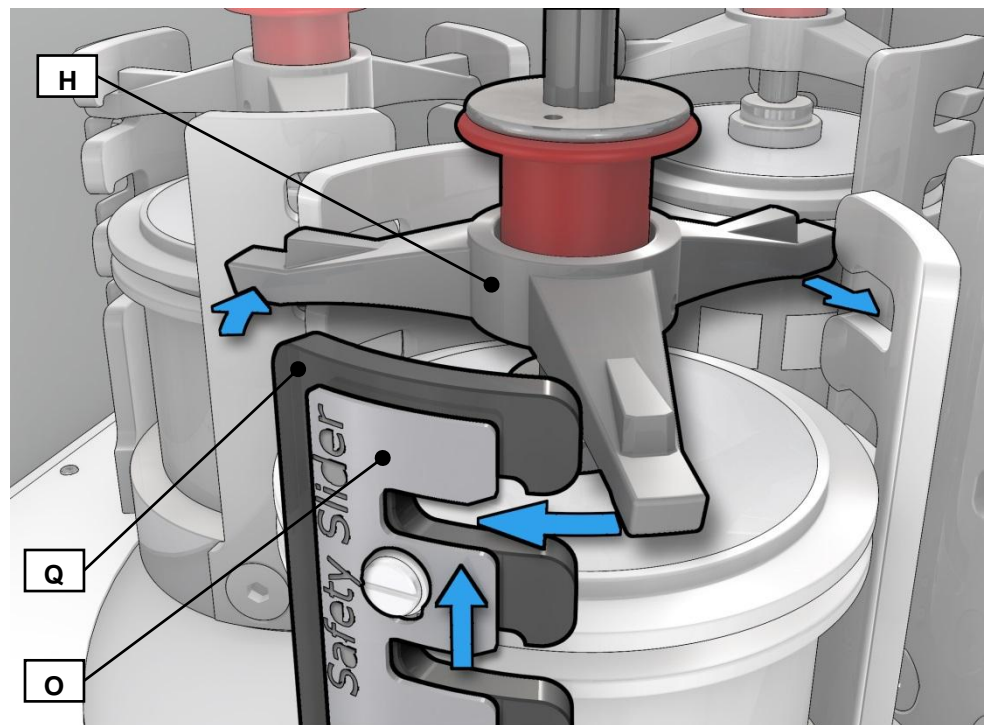


Fig. 19: Inserting the clamping unit

- Insert the clamping unit (H) in the three latching brackets (Q).

5.8.4 Function of the locking sleeve

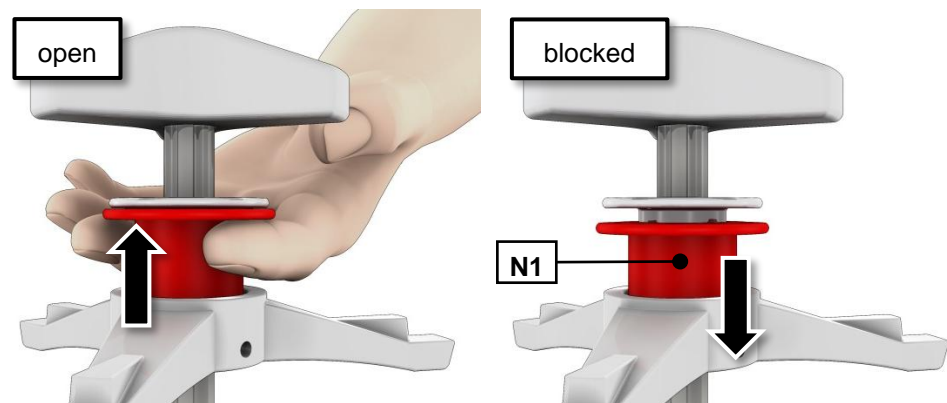


Fig. 20: Function of the locking sleeve

- Pull the red sleeve (N1) upwards and clamp the grinding jar by twisting the three-star grip to the right.
- Allow the red sleeve (N1) to audibly click down into place, if necessary tightening slightly using the three-star grip.
- It should no longer be possible to turn the three-star grip.
- The locking sleeve now in the blocked position stops the threaded spindle from coming loose automatically.

5.9 Releasing the grinding jar clamping mechanism

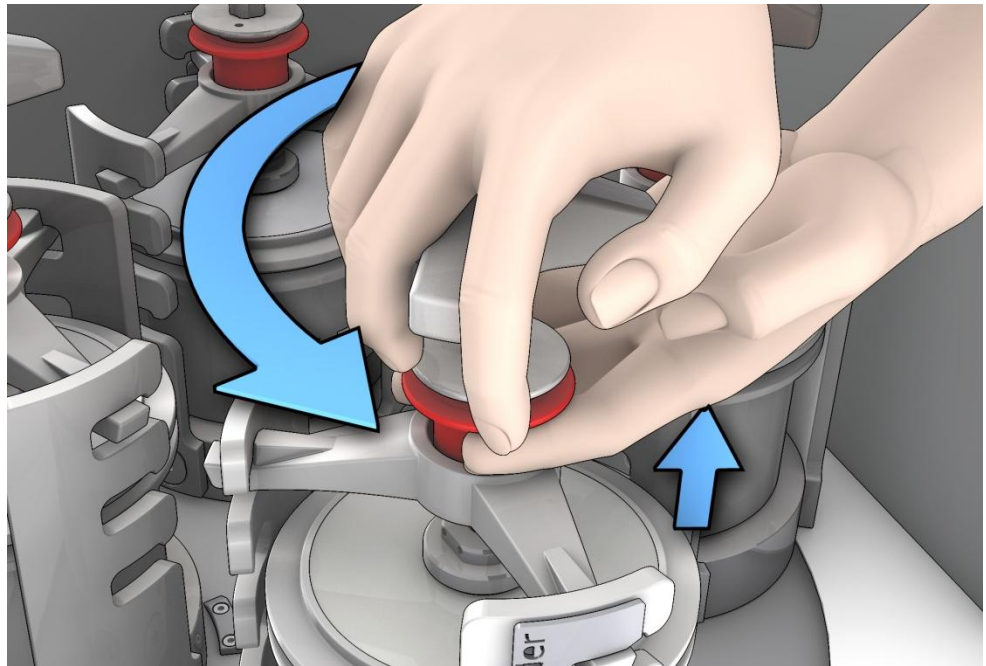


Fig. 21: Releasing the grinding jar clamping mechanism

- Pull the red sleeve (5) upwards.
- Turn the three-star grip to the left to release the grinding jar.
- Keep turning the three-star grip to the left until the grinding jar clamping mechanism can be removed.

5.10 Opening the Clamping Mechanism with the Opening Aid for the Clamping Unit

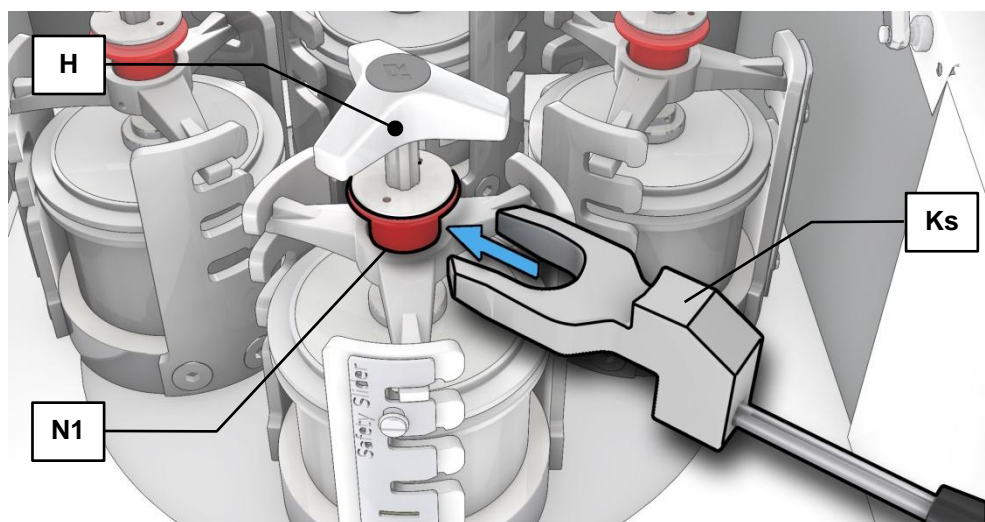


Fig. 22: Inserting the opening aid for the clamping unit

- The opening aid for the clamping unit (**KS**) [accessory] can be used to loosen or tighten the clamping unit (**H**).
 - When inserting the opening aid for the clamping unit (**KS**), the locking sleeve (**N1**) is pushed upwards and released.

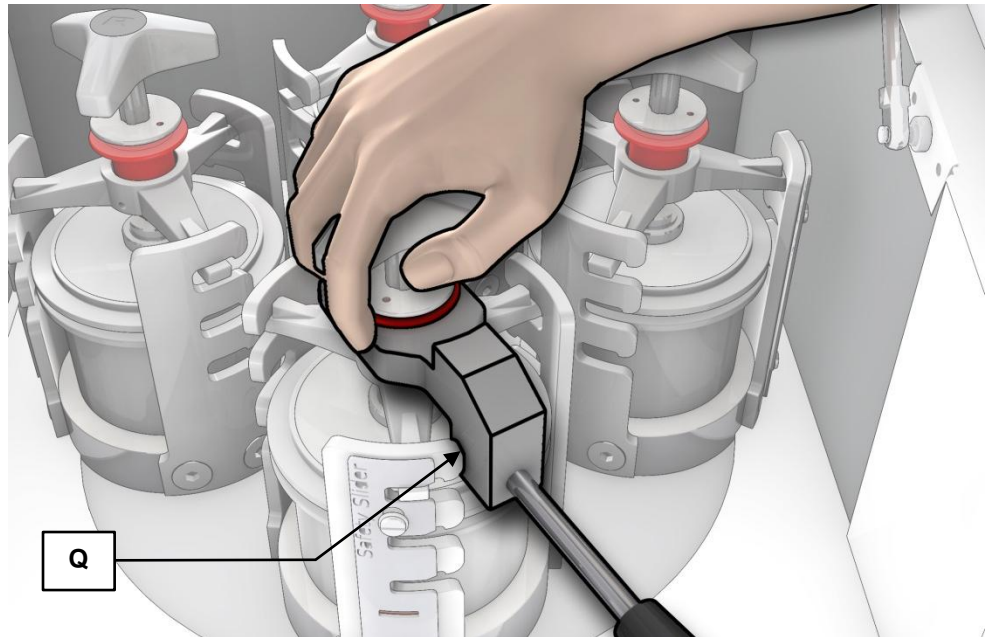


Fig. 23: Loosening the clamping unit

- Rest the opening aid for the clamping unit (**KS**) on the latching bracket (**Q**) during closing and loosening.

5.11 Power failure during grinding

If the power fails during grinding, grinding will be interrupted. All parameters are retained, and the remaining grinding time is saved. When the device is switched on again, a corresponding message appears in the display.

You can continue the process by pressing the START button. The automatic saving of the remaining grinding time means the grinding then continues until the end of the originally set grinding time.

For safety reasons the device must be opened and the grinding chamber inspected. Grinding then continues automatically when the lid is closed.

The process is aborted by pressing the STOP button.

5.12 Mahlbehälterauswahl für unterschiedlichen Probenmaterialien

This device is only suitable for use with grinding jars from Retsch GmbH with a nominal volume of 12 ml - 500 ml.

They are available in the following materials:

- Agate
- Sintered corundum
- Zirconium oxide
- Stainless steel
- Special steel
- Tungsten carbide

5.13 Sample quantity

NOTICE

If the fill level in the grinding jar is too high or too low, this will impair the milling result and may cause damage (increased abrasion) to the grinding set.

5.13.1 Guides for material quantity and balls

Grinding jar volume	Sample volume	Max. feed size	Ball filling (each)					
			Ø 5 mm	Ø 7 mm	Ø 10 mm	Ø 15 mm	Ø 20 mm	Ø 30 mm
12 ml	≤ 5 ml	< 1 mm	50	15	5	-	-	-
25 ml	≤ 10 ml	< 1 mm	95-100	25-30	10	-	-	-
50 ml	5 - 20 ml	< 3 mm	200	50-70	20	7	3-4	-
80 ml	10 - 35 ml	< 4 mm	250-330	70-120	30-40	12	5	-
125 ml	15 - 50 ml	< 4 mm	500	110-180	50-60	18	7	-
250 ml	25 - 120 ml	< 6 mm	1100-1200	220-350	100-120	35-45	15	5
500 ml	75 - 220 ml	< 10 mm	2000	440-700	200-230	70	25	8

In addition to the instrument settings, the filling level of the grinding jar is also of crucial importance for a successful grinding process in the Planetary Ball Mill of Retsch GmbH. The usable capacity of the grinding bowls depends on the type of material. The specified number of balls is the minimum quantity per grinding bowl. An optimal grinding result is achieved with the larger number of permissible grinding balls, if specified. In exceptional cases, the number of grinding balls can be reduced by up to 15 %, but then increased abrasion of the grinding set must be expected.

When grinding bulk materials, the grinding jar filling should consist of approximately one third of sample and one third of ball quantity. The remaining third is the free grinding jar volume, which is required for the movement of the balls.

If an increase or decrease in sample volume is to be expected during the grinding process, the amount of sample can be adjusted within the range listed in the table. Thus, e.g. for voluminous materials such as wool, leaves, grasses and similar, a material filling level of 70 – 80 % is necessary. For wet grinding with grinding balls < 3 mm, the ball charge should be 60 % of the grinding jar volume.

5.14 Ultrafine grinding

High degrees of fineness can frequently only be achieved using wet grinding.

Improved final fineness can be achieved in dry grinding by adding a few drops of stearin or acetic acid and by using grinding balls with a $\text{Ø} < 10 \text{ mm}$ and a fill level of 70-80% of the grinding jar volume.

5.15 Stacking the grinding jars

Stacking of grinding jars is possible in the sizes 12 ml, 25 ml, 50 ml and 80 ml.

A maximum of two grinding jars may be stacked on top of each other.

In order to insert and stack specific grinding jars, you may need the adapters which are available as accessories.

5.15.1 Stacking the 50 ml and 80 ml grinding jars

50 ml grinding jars in Comfort design and 50 ml and 80 ml grinding jars in EasyFit design can be stacked. You may need the adapters which are available as accessories for stacking.

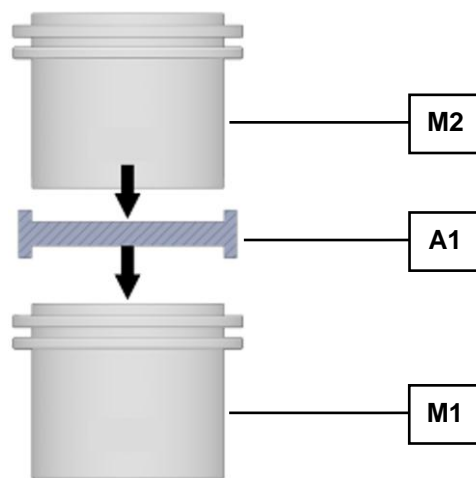


Fig. 24: Stacking grinding jars

Stacking process:

- Place the grinding jar (**M1**) in the grinding jar plate
- Place the adapter (**A1**) on the lid of M1
- Place grinding jar (**M2**) on the adapter
- Clamp the grinding jars as described in the chapter “Clamping the grinding jars”.

5.15.2 Stacking the 25 ml and 12 ml grinding jars

Grinding jars with a capacity of less than 50 ml can be stacked on top of each other directly without the use of accessories.

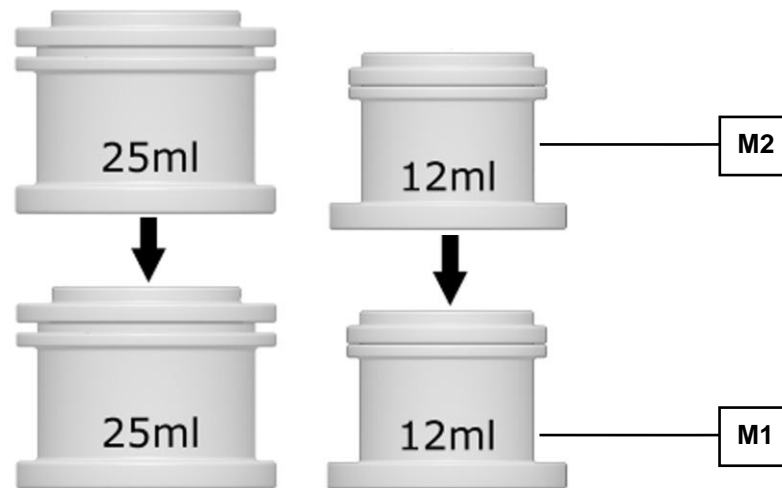


Fig. 25: Stacking the 25 ml and 12 ml grinding jars

Stacking process:

- Place the grinding jar (**M2**) directly on top of grinding jar (**M1**).
- Clamp the grinding jars as described in the chapter “Clamping the grinding jars”.

5.16 Handling grinding jars

5.16.1 Carrying and gripping



Fig. 26: Grinding jar

The gripping flanges (**Gr**) on the grinding jar lid permit secure handling.


5.16.2 Anti-rotation device

All 250ml and 500ml grinding jars have a drilled hole underneath the grinding jar as anti-rotation device. The anti-rotation device is only used with the PM400 and is located on the side of the grinding jar casing with lettering.

5.16.3 Heating the grinding jars

Depending on the grinding time and the fill level, the grinding jars can be heated to a temperature of up to 150°C during grinding.

This temperature change causes the pressure inside the grinding jar to increase. When unscrewing the lid, please note that this excess pressure is released by a sudden escape of air. This may be accompanied by particles of sample material.

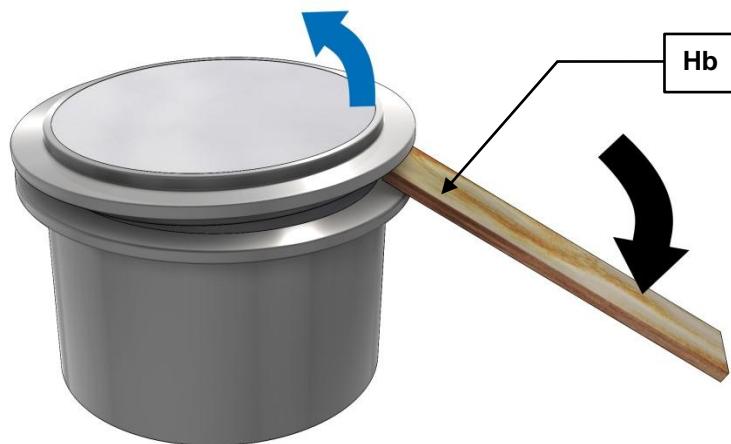

When grinding jars have been left to cool, a vacuum is created inside them which can make them difficult to open. The grinding jars can be prised open between the gripping flanges on the lid and the grinding jar, e.g. using a flat wooden stick (**Hb**).

5.17 Grinding jar identification

All grinding jars can be identified by lettering which shows the article number and material.

5.18 Grinding jar cleaning

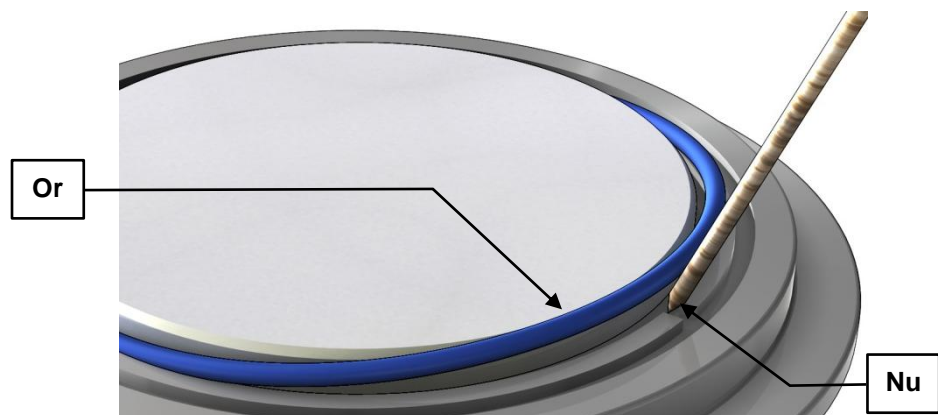


Fig. 27: Removing the O-ring

The O-ring (**Or**) on the groove (**Nu**) on the underneath of the lid can be lifted out easily to clean the grinding jar.

Grinding jars, including those with glued ceramic inserts, can be cleaned using alcohol, benzine or normal household detergents

NOTICE

When cleaning, do not expose grinding jars with ceramic inserts to sudden temperature differences.

Sudden temperature fluctuations can cause the ceramic inserts to crack.

5.18.1 Drying the grinding jars

After cleaning, the grinding jars can be dried in the drier cabinet at the temperatures specified below.

Material of the grinding jars	Temperature
Hardened stainless steel	up to 200°C
Tungsten carbide	up to 120°C
Sintered corundum	up to 120°C
Agate	up to 120°C
Zirconium oxide	up to 120°C

5.19 Opening and closing of the grinding jars with safety locking device

A safety locking device is recommended for grinding processes (especially wet grinding) where an increase in pressure is to be expected in the grinding jar. If a gassing lid is used, the grinding jar must also be sealed with a safety sealing device.

NOTE

For Comfort design grinding jars, the safety sealing device must always be selected taking into account the grinding jar size and material.

The use of agate grinding jars for wet grinding with solvents should be given particular consideration due to the internal pressures that arise and the non-homogeneous material properties of a natural product.

The type of safety locking device must be adapted to the grinding jar used. The safety locking device with two clamping rings must be used for grinding bowls in the Comfort design and with locking hooks for grinding bowls in the EasyFit design. No product combinations other than those described here may be used!



Fig. 28: A Comfort design grinding jar (left) and an EasyFit design (right)



Fig. 29: Safety locking device for Comfort design grinding jars

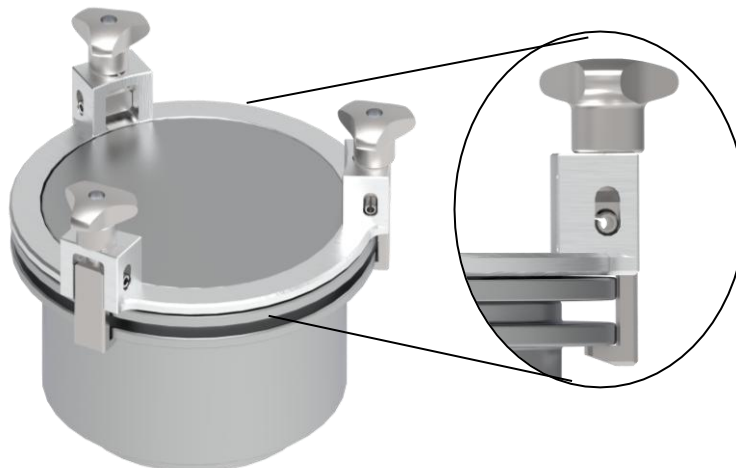


Fig. 30: Safety locking device for EasyFit design grinding jars

Tighten the screw connection on the safety locking device alternately and gradually. Tightening unevenly can cause the cover and the safety locking device to tilt!

With the safety locking device for grinding jars in the Comfort design, the three clamping screws must be tightened to 2.5 Nm.

With the safety locking device for grinding jars in the EasyFit design, the three clamping screws (using the opening and closing aid) must be tightened to at least 8 Nm.

Internal pressures up to a max. 5 bar are only permissible with this pre-tensioning.

NOTE

Damage to the grinding jar lid and the appliance.

The three screws on the safety clip of the gassing cover can loosen and damage the inside of the appliance.

After tightening the grinding jar clamping device, check whether the three screws on the safety locking device are still tight.

NOTE

The appliance is equipped with a fan that extracts the waste heat generated during grinding directly from the grinding chamber. The extraction volume per hour is more than 20 times the grinding chamber volume. The fan has a standstill monitor with signalling.

If necessary, divert the air flow from the fan into a fume cabinet during grinding.

Before removing the grinding jar, check that the safety locking device is firmly in place.

Only remove the grinding jar with the safety locking device and only open it in a safe position (suction device) after it has cooled down.

5.20 Wet Grinding with Highly Flammable Materials

Wet grindings using highly flammable materials are permissible with this device when complying with specific precautionary measures.

When using highly flammable materials as grinding aid, such as hexane, isopropanol, ethanol, petrol or similar, it is to be assumed that the inside of the grinding jars should be classified as zone 0, i.e. a constantly present explosive mix.

Therefore, it is necessary to prevent potentially explosive vapours during the grinding process from escaping from the clamped grinding jars and from entering areas in which the requisite

ignition energy is present. These vapours may be pressed out of the grinding jar in particular by the heating which takes place and the resulting pressure increase inside the grinding jar.

It is therefore urgently recommended that the managing operator (employer) using the device assesses the dangers that exist before using corresponding solvents, taking local conditions into account, as part of a coherent explosion protection concept. If necessary, supplementary organisational measures should be recorded in writing in an explosion protection document.

This procedure is regulated in the EU under Articles 118 and 118a of Directive 89/391/EEC. In other countries outside the EU, the comparable provisions are to be observed.

The following must be verified with respect to the device:

- **Only grinding jars with safety closure devices may be used!**
- It is necessary to consider the durability of the O-rings (EPDM 75° shore) when selecting the solvent and the durability of the adhesive used when using ceramic inserts.
- The safety closure devices of the grinding jars must be tightened firmly.
- Please note that the grinding jars can heat up considerably depending on the grinding jar size, the ball filling, the speed and the grinding duration.
- Before removing the grinding jars, the tightness of the safety closure devices must be checked anew.

6 Display and operation

6.1 Symbols in the Display Unit

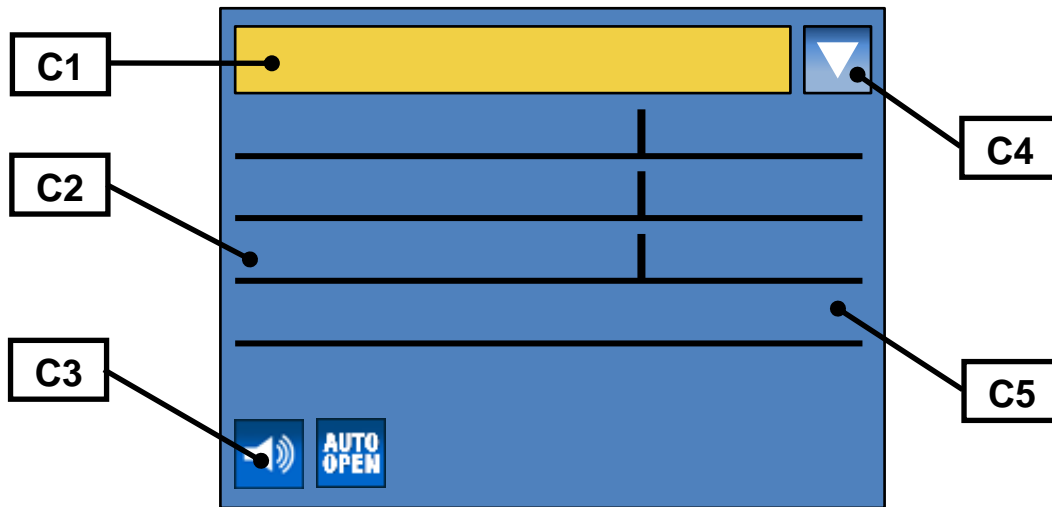


Fig. 31: View of the menu on the display unit

Element	Description	Function
C1	Menu navigation	Switching between manual operating mode, program and basic settings
C2	Specification of grinding parameters	Displaying and setting grinding parameters
C3	Icons for device functions	Displaying the function statuses of sound, automatic opening and grinding jar recognition
C4	Icon for scrolling direction	Displays the possible scrolling directions
C5	Grinding parameters	Display of values

	Automatic opening switched on
	Automatic opening switched off
	Direction reversal switched on
	Direction reversal switched off
°C	Motor or frequency converter too hot
Power	Grinding output display
	Acoustic warning signal on
	Acoustic warning signal off
	Scrolling upwards or downwards possible
	Only scrolling upwards possible
	Only scrolling downwards possible

6.2 Display unit – operation of the device

This device offers a new, very convenient user interface. All relevant data can be entered and retrieved using a graphics display with one-button operation. The menu is available in different languages.

6.2.1 Adjustment options using the display menu

The selection bar in the display should be operated as follows:

Rotating function I)

- Rotate the operating button to get to the different menu items. The selected menu items are marked by the dark selection bar. Areas that cannot be changed are skipped.

Rotating function II)

- Rotate the operating button to change numerical values and decisions in the menu items.

Press I)

- Press the operating button to open selected menu items.

Press II)


- Press the operating button to confirm settings.

Press III)

- Keeping the operating button pressed for longer takes you back to the basic screen (Level 1). Navigating between operating modes
- Rotate the operating button in a clockwise direction until the dark line cursor is in the navigation menu (**C1**).
- Press the operating button (**F**).
- The icon for the scrolling direction (**C4**) changes from




- By pressing the operating button, navigate between manual operation, Programs 01 to 10 and the basic settings operating modes.
- Press the operating button (**F**) to activate the selected operating mode.
- The icon for the scrolling direction (**C4**) changes from

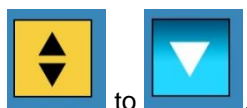

- By rotating the operating button, switch to the sub-items of the selected menu item.

6.3 Direct access to the language menu

If you have unintentionally set the wrong language, you can go straight to the language menu by following the steps below.

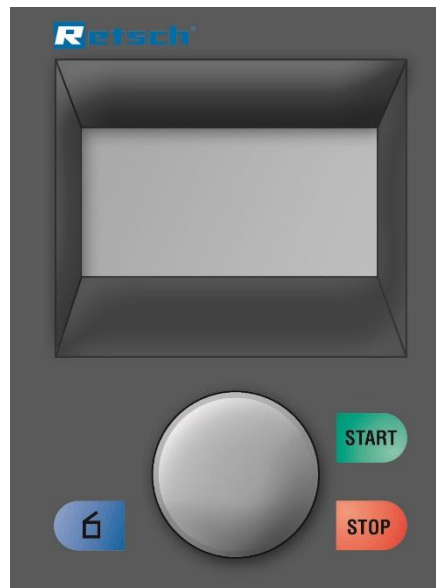


- Switch the device off at the main switch.
- Switch the device on while simultaneously pressing the buttons **START - STOP – Open hood**.
- After selecting the correct language, switch the device off and then immediately back on.
- Confirm your selection by pressing the operating button.

The device is now set permanently to your language and you are in the main menu.

6.4 Menu structure

Complete summary of all menu options:

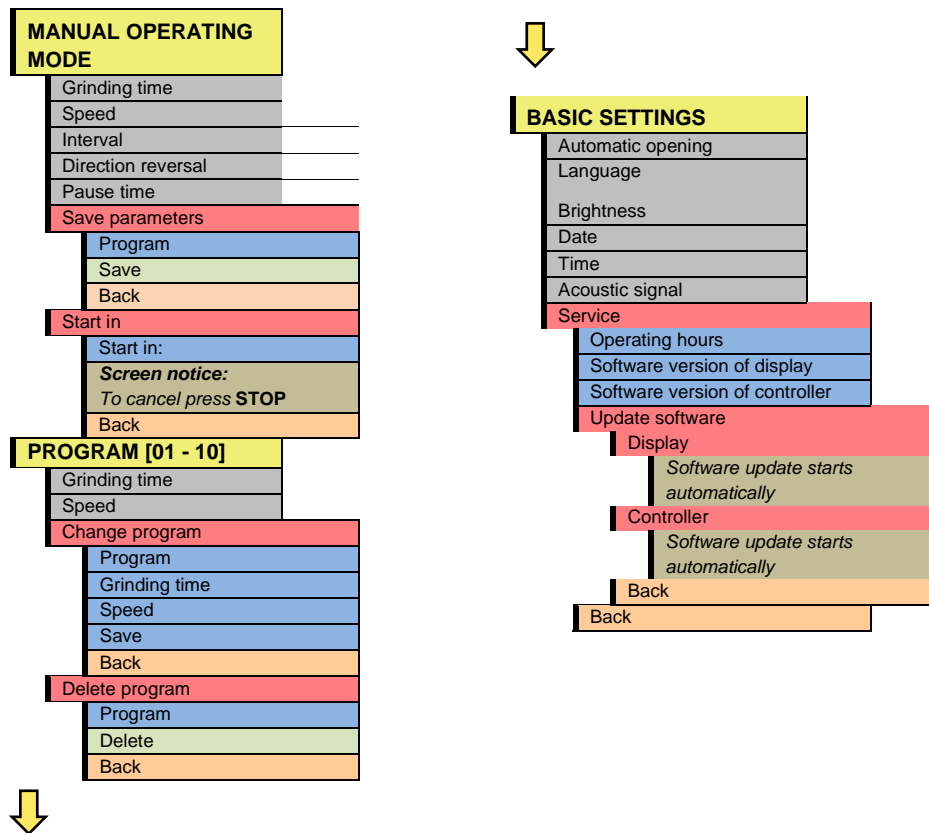

6.5 Operating modes

You can select the following operating modes using the menu navigation (**C1**):

6.5.1 Manual operation

When this function has been set you can access and change all parameters and functions at any time. This is also possible during grinding.

6.5.2 Program 01 to 10

In programs 01 to 10 the previously set parameters for grinding time and speed can be saved to a memory.

6.5.3 Basic settings

You can adjust the following device settings in this settings menu:

- Automatic opening
- Language
- Brightness
- Date
- Time
- Acoustic warning signal
- Service

6.6 Manual Mode

6.6.1 Grinding time

00:00:01 to 99:59:59 (hours : minutes : seconds)

The device is started with the preselected grinding time and at the last speed used. Direction reversal with pause time has not been switched on

6.6.2 Speed

30 to 400 rpm

The device is started with the preselected grinding time and at the last speed used. Direction reversal with pause time has not been switched on

6.6.3 Interval

00:00:01 to 99:59:59 (hours : minutes : seconds)

The interval time can be set here according to the grinding time. If no interval has been set, no direction reversal can be set.

6.6.4 Direction reversal

On/off

The device is started with the preselected grinding time, speed and direction reversal. The machine rotates in one direction for the set interval time, comes to a halt and starts again immediately in the other direction without a pause time.

6.6.5 Pause time

00:00:01 to 99:59:59 (hours : minutes : seconds)

The pause time between intervals can be set here.

No pause time can be set if no interval has been set.

The device is started with the preselected grinding time, speed and direction reversal. The machine rotates in one direction for the set interval time and comes to a halt; once it has stopped at the set interval time, the previously set pause time is displayed and counts down to 00:00:00.

The device starts again in the other direction when the pause time has expired.

6.6.6 Save parameters

The previously set parameters such as grinding time and speed can be saved in a memory here.

- Set the desired parameters.
- By turning the control knob (F), switch to the **Save parameters** menu option.
- Press the control knob (F).
- The **Save parameters** menu opens and the dark cursor line is on **Program**.
- Press the control knob (F) to select a program memory location.
- By turning the control knob (F), switch to the desired memory location.
- Press the control knob (F) to exit the memory location selection.
- Select either
 - **Save** to save settings or
 - **Back** to cancel without saving.

6.6.7 Start in

00:00:01 to 99:59:59 (hours : minutes : seconds)

You can set a countdown for starting the device.

- Press the STOP button to cancel the countdown.

6.7 Programme Mode

6.7.1 Change program

You can change the saved parameters for any program in this menu.

- By turning the control knob (F), switch to the **Change program** menu option.
- Press the control knob (F).
- The **Save parameters** menu opens and the dark cursor line is on **Program**.

NOTICE

You can change the active program or any other program.

- Press the control knob (F) to activate the program selection.
- By turning the control knob (F), switch to the desired memory location.
- Press the control knob (F) to exit the memory location selection.
- Adjust the desired grinding parameters.
- Finally select either
 - **Save** to save settings or
 - **Back** to cancel without saving.
- This takes you back to the Program level.

NOTICE

It is not possible to start a program that has not been saved.

6.7.2 Delete program

You can delete the saved parameters for any program in this menu

NOTICE

Only the saved parameters for the program concerned are deleted. The program memory location continues to exist.

- By turning the control knob (F), switch to the **Delete program** menu option.
- Press the control knob (F).
- The **Delete program** menu opens and the dark cursor line is on Program.
- Press the control knob (F) to activate the program selection.
- By turning the control knob (F), switch to the desired program.
- Press the control knob (F) to exit the program selection.
- Finally select either
 - **Delete** to delete the settings or
 - **Back** to cancel without deleting.
- This takes you back to the Program level.

6.8 Basic settings

NOTICE

It is not possible to begin grinding while the basic settings menu is active.

6.8.1 Automatic opening

In this menu you can set whether the grinding jar lid opens automatically when grinding ends or only when the button is pressed.

If the function is switched off, the following pictogram appears on the display as confirmation.



Fig. 32: Automatic opening pictogram

6.8.2 Language

You can select the menu language here. After selecting and pressing the control knob, the entire menu structure is displayed in your language.

NOTICE

The language menu is displayed when the device is switched on for the first time.

- Select the desired language by turning the control knob.
- Press to confirm the selection; “open lid” appears on the display.

6.8.3 Brightness

The brightness can be adjusted to suit the respective user or environment (sunshine, glare etc.).

6.8.4 Date

The current date can be entered here.

The device can be disconnected from the mains for up to 30 days before the settings are lost.

6.8.5 Time

The time can be entered here.

The time then appears in the stand-by monitor.

The device can be disconnected from the mains for up to 30 days before the settings are lost.

6.8.6 Acoustic warning signal

Error messages indicating incorrect operation can be supported by an acoustic warning signal.

The corresponding pictogram appears if the function has been switched off

6.8.7 Service

6.8.7.1 Operating hours

Grinding hours are counted, i.e. the total times between START and STOP. It is not possible to manipulate the times.

6.8.7.2 Software version of display

Shows the software version of the display.

6.8.7.3 Software version of the controller

Shows the version of the operating software .

6.8.7.4 Update software

SOFTWARE VERSION

The version of the operating software can be called up and updated as required. If necessary, get in touch with your Retsch distributor.

Should you have reached this menu by mistake and cannot return to the previous menu, switch the device off at the main switch and re-start it.

6.8.7.4.1 Safety notice

For many years the grinding jar clamping mechanism has been a proven, easy to use and reliable device. The basic prerequisite both for the safety of the operator and for the life of the machine components is careful clamping of the grinding jars.

Please remember that this device involves grinding equipment which applies very high amounts of energy into the sample material, and that the grinding set therefore needs to be locked into position carefully.

Correct securing of the grinding jar is always queried prior to starting the machine in order to prevent operating errors.

Employees who have been especially trained and who are familiar with operation of the PM can also permanently fade out this safety notice. We do not recommend this course of action, particularly when there are changes in operating staff!

The device software is set up such that the message about the need to confirm clamping of the grinding jars is always displayed when the start button is actuated prior to starting the machine.

The grinding process begins after confirmation.

This safety notice can be faded out in the "Settings" menu.

7 Fault messages

Error code	DEFECT DESCRIPTION TRANSLATION
E10	DRIVE OVERLOAD
E20	FAILURE MAIN BOARD
E23	FAILURE FAN
E25	FAILURE DISPLAY
E26	FAILURE FREQUENCY CONVERTER
E41	FAILURE SPEED SENSOR
E50	FAILURE IN SAFETY CIRCUIT
H10	ALLOW DRIVE TO COOL DOWN
H13	OVERLOAD! REDUCE SPEED!
H14	OVERLOAD! SPEED HAS BEEN REDUCED!
H42	OPEN AND CLOSE LID/COVER
H45	MAINS INTERRUPTION

8 Installing additional equipment

By means of the Retsch additional equipment, the planetary ball mills PM 100, PM 300 and PM 400 can be flexibly adapted to different working conditions.

The following additional equipment is available for the Retsch planetary ball mills PM 100, PM 300 and PM 400:

- Adapter for 24 x 1.5 ml glass vials
- Adapter for 7 x 20 ml glass vials

⚠ CAUTION If the additional equipment is loaded incorrectly, the grinding vessels may splinter during the grinding process. Please observe the specifications regarding the sizes of the grinding balls and the maximum permitted speeds.

8.1 Adapter for glass vials

NOTICE

Strong vibration and loud noise

Uneven load

- The device can produce very high levels of vibrations and noise when loading is uneven.
- **Always use 2 grinding jars opposite each other.**
- **The grinding stations must be operated with the identical grinding jars and with the same weight during each grinding process.**
- **Switch the device off immediately if there are high levels of vibration and noise and check the number and gross weight of the jars.**

The device can be equipped with an adapter for glass vials. The adapters enable the simultaneous use of up to 24 x 1.5 ml or 7 x 20 ml samples in the respective vials.

Mount the adapter as follows:

NOTICE Always fit the adapter in a symmetrical arrangement. Otherwise, the grinding jar lid may tilt and the adapter may not be clamped properly.

⇒ Insert the compressure spring into the desired opening of the adapter.

⇒ Place the loaded and closed glass vial on the compressure spring with the lid facing upwards. The tension of the spring pushes the vial upwards.

⇒ Insert the loaded adapter into the grinding jar support of the desired device.

Close the adapter with the lid and secure it with the clamping unit. It must be ensured that the lid sits straight and is firmly clamped in place.

⚠ WARNING Please observe the instructions in the operating instructions for your device regarding the safe handling of the clamping unit. The lid must close the adapter flush and must not rest at an angle!

⇒ Alternatively, the lid of the adapter can also be secured with the safety locking device for grinding jars. In such a case, load the adapter as described above, put on the lid and tighten the safety locking device. Subsequently, insert the closed adapter into the grinding jar support.

NOTICE In order to reduce wear and tear on the compressure springs, it is recommended that only springs are inserted into the openings of the adapter which are loaded with vials.

⚠ CAUTION The clamping force of the compressure springs pushes the inserted vials out of the adapter. Therefore, transport the adapter which is loaded with vials without the lid. Alternatively, secure the lid with the safety locking device.



Fig. 33: Exploded view of the adapter with compressure spring and 1.5 ml vial



Fig. 28: Permissible arrangement of the glass vials in the adapter

⚠ CAUTION Glass vials are not suitable for grinding hard-brittle materials. These are disposable glass vials which must not be used more than once. Please observe the maximum speed specifications.

⚠ CAUTION Glass vials can break and cause cuts!

The filling of the vials deviates specifically from the 1/3 rule in terms of application and filling:

Example 1 (1.5 ml)	
Grinding balls	2 – 4 x 3 mm grinding balls, stainless steel
Sample material/ Mixture of reactants	20 % filling volume of the vial, top up with a few µl of solvent

Example 2 (1.5 ml)	
Grinding balls	approx. 1/3 of the glass vial with 3 mm grinding balls, stainless steel
Sample material/ Mixture of reactants	max. 20 % filling volume of the vial, top up with a few μ l of solvent

Example 3 (20 ml)	
Grinding balls	approx. 1/4 of the glass vial with 3 mm grinding balls (stainless steel, zirconium oxide, agate)
Sample material/ Mixture of reactants	50 % filling volume of the vial, top up with a few μ l of solvent

8.1.1 Speed limits

Due to the high energy input of the device, strong forces act on the glass vials. The maximum speed specifications must be adhered to in order to prevent any damage to the adapter and the vials.

Device	Maximum speed	
	Adapter for 24 x 1.5 ml glass vials	Adapter for 7 x 20 ml glass vials
PM 100	550 rpm	350 rpm
PM 300	500 rpm	300 rpm
PM 400	400 rpm	250 rpm

8.1.2 Energy input

The energy input of the device to the glass vials in the adapter differs depending on the arrangement of the openings on the inner or outer diameter. The force applied to the inner diameter is comparable with 50 - 125 ml grinding jars. The force applied to the outer diameter is comparable with 250 - 500 ml grinding jars.

Green: Inner diameter, 50 - 125 ml grinding jars

Red: Outer diameter, 250 - 500 ml grinding jars

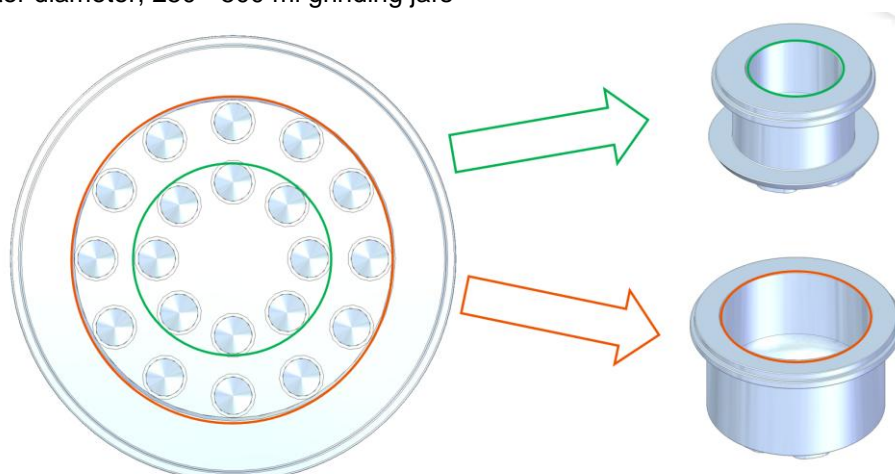


Fig. 34: Energy input as a function of diameter

9 Cleaning, Wear and Maintenance

WARNING

Risk of a fatal electric shock

- An electric shock can cause injuries in the form of burns and cardiac arrhythmia, respiratory arrest or cardiac arrest.
- **Do not clean the blender under running water. Use only a cloth dampened with water.**
- **Disconnect the power supply plug before cleaning the blender.**

WARNING

The device must always be switched off and disconnected from the mains before any interventions for cleaning or servicing purposes.

W0012

9.1 Maintenance

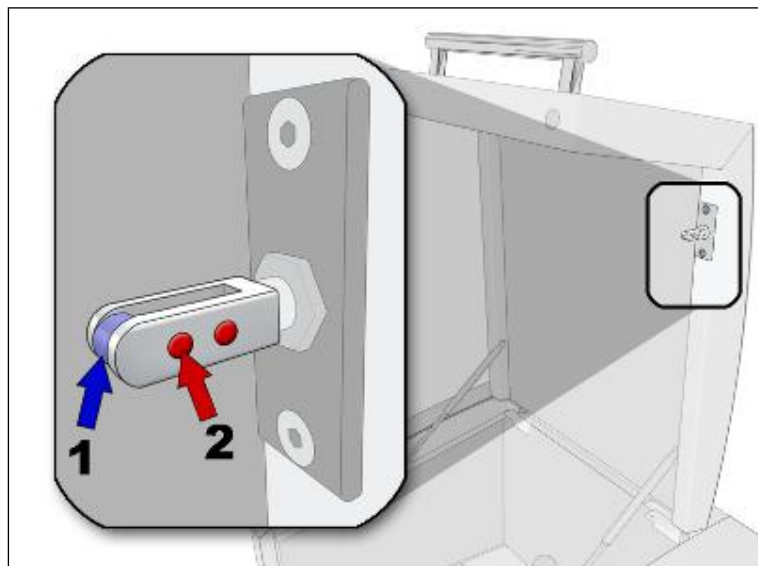


Fig. 35: Closing pin maintenance

The following maintenance work should be carried out at regular intervals, and at least monthly to guarantee the operational safety of the device:

- Check roller (1) of the closing pin for free movement, and oil where necessary, e.g. using sewing machine oil.
- Clean the magnets (2) on the closing pin.

9.1.1 Servicing the clamping unit

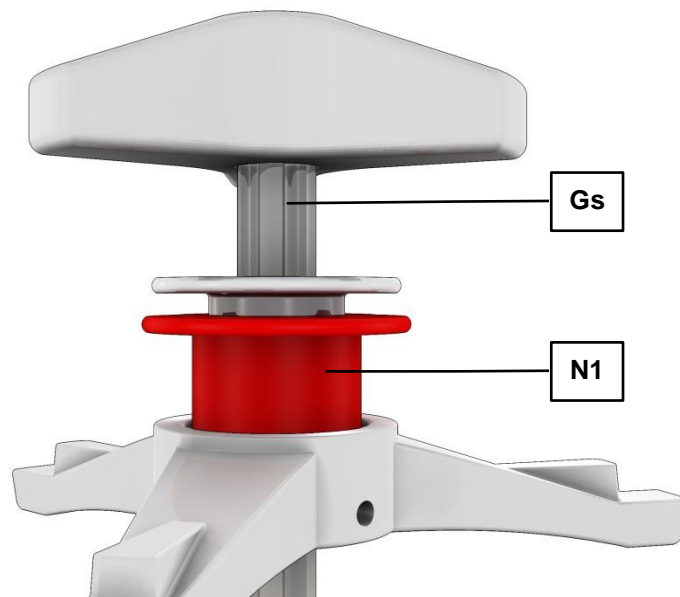


Fig. 36: Servicing the clamping unit

- Check that the threaded spindle (**Gs**) and locking sleeve (**N1**) can move freely.
 - In most cases a drop of oil will help.

The smooth running of the threaded spindle and locking sleeve is the prerequisite for the secure clamping of the grinding jars.

Locking sleeves which do not slide down automatically through spring force cannot reliably prevent the threaded spindle from coming loose. The grinding jars may be ejected.

9.1.2 Wear to the bases of the clamping unit

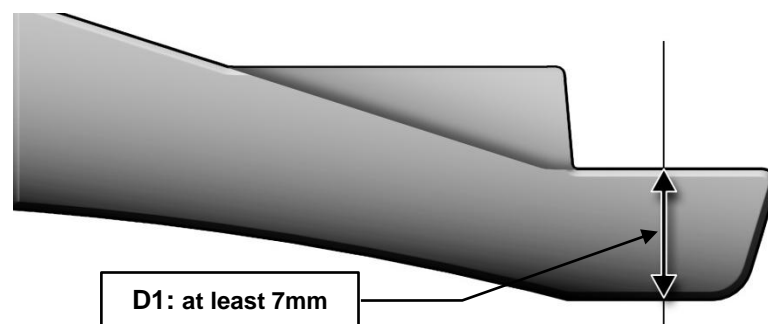


Fig. 37: Wear to the bases of the clamping unit

- Check the thickness (**D1**) of the three lead frame bases regularly for wear (at least monthly).
 - The thickness (**D1**) of the three lead frame bases must not be less than 7mm.
 - Operational safety can no longer be guaranteed if the thickness falls below this value (**D1**). The grinding jars may be ejected.

9.1.3 Rubber washer on the pressure plate

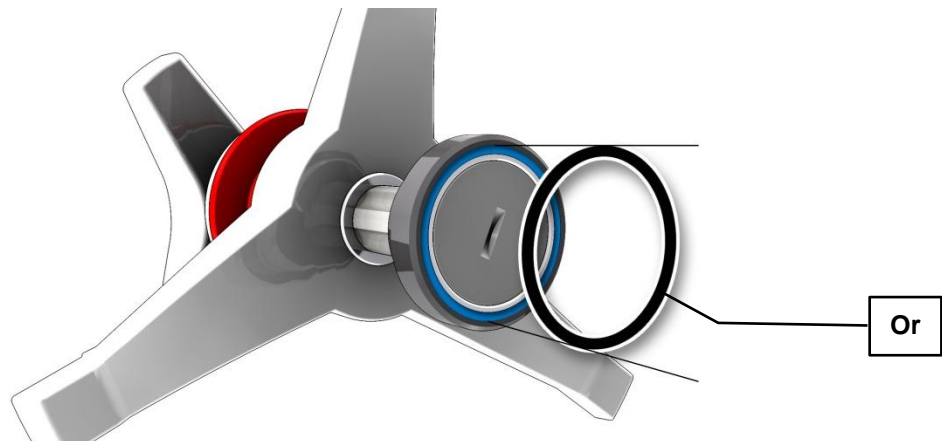


Fig. 38: Rubber washer on the pressure plate

- Check the O-ring (**Or**) on the pressure plate regularly for wear and to ensure it is secure.

9.1.4 Wear to latching bracket

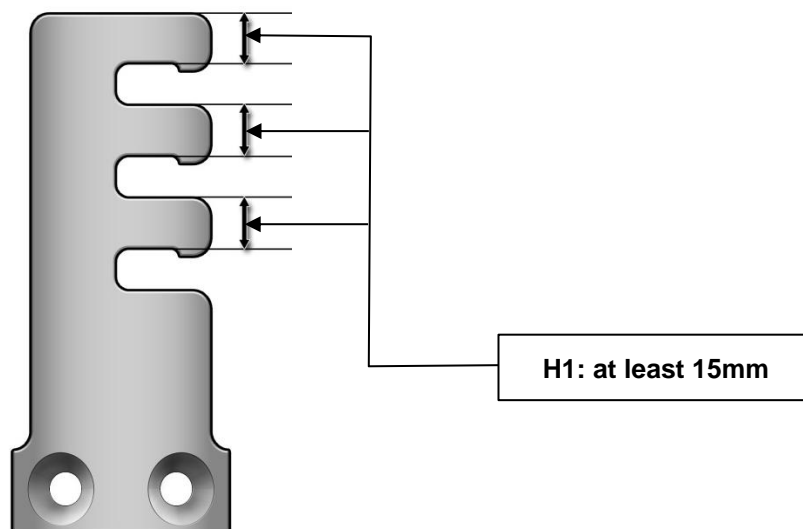


Fig. 39: Wear to the latching bracket

- Check the thickness (**H1**) of the twelve latching brackets (**Q**) regularly for wear (at least monthly).
 - The thickness (**H1**) of the twelve latching brackets (**Q**) must not be less than 15mm.
 - Operational safety can no longer be guaranteed if the thickness falls below this value (**H1**). The grinding jars may be ejected.

10 Return for Service and Maintenance



Fig. 40: Returned goods dispatch note

RETSCH devices and accessories can only be accepted for repair, maintenance or calibration if the returned goods despatch note has been correctly completed in full.

- When returning a device, attach the returned goods dispatch note to the outside of the packaging.

In order to eliminate any health risk to our employees, we reserve the right to refuse acceptance and to return the respective delivery at the expense of the sender.

11 Disposal

Please observe the respective statutory requirements with respect to disposal.

Information on disposal of electrical and electronic machines in the European Community.

Within the European Community the disposal of electrically operated devices is regulated by national provisions that are based on the EU Directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE).

Accordingly, all machines supplied after 13.08.2005 in the business-to-business area to which this product is classified, may no longer be disposed of with municipal or household waste. To document this they have the following label:

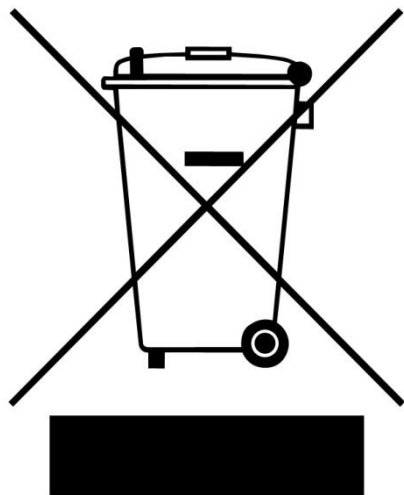


Fig. 41: Disposal label

Since the disposal regulations within the EU may differ from country to country we would request you to consult your supplier.

12 Index

1	
16A.....	20
A	
Acoustic warning signal	49
Adapter	51
adapter for 24 x 1.5 ml glass vials	51
adapter for 7 x 20 ml glass vials	51
Adjustment options using the display menu	43
Agate.....	34
all current sensitive	14
all current sensitive	14
Ambient temperature	12
Amperage	12
Anti-rotation device	37
Automatic lid closing	21
Automatic opening	48
B	
Ball	
quantity.....	35
size	35
Ball filling.....	35
Bar code	12
Basic settings.....	46, 48
Brightness	48
C	
Capacity	12
CE marking	12
Ceramic insert.....	41
Change program	47
Changes	6
Clamping bolt.....	28
Clamping mechanism	
opening	33
Clamping unit	
servicing.....	55
wear.....	55
Cleaning.....	54
Cleaning the grinding jars	38
Clearance distance	22
Close-up of the grinding chamber	23
Closing pin	25
Confirmation	10
Connection cable	14, 21
Copyright	6
creating interface connection.....	14
D	
Date	48
Decrease of sample volume during grinding	
process.....	35
Degree of protection	21
Delete program	48
Depth	22
Description.....	24, 25, 42
Device	
closing	25
open.....	25
Device designation.....	12
Dimensions	22
DIN 45635-31-01-KL3.....	21
Direct access to the language menu	44
Direction reversal	46
Display and operation	42
Display unit – operation of the device.....	43
Disposal	58
Disposal label.....	12, 58
E	
Electrical connection	14, 20
EMC Directive	14
Emergency release	26
Emergency unlocking.....	26
Emissions.....	21
Energy input.....	53
Error code	50
Error messages.....	50
Errors	50
Explanations of the safety instructions	7
External fuse	14, 20
F	
Fault messages.....	50
Feed grain size	20
Feed sizes.....	35
Frequency converter.....	14
Function	24, 25, 42
Function of the locking sleeve	32
Fuse strength	12
Fuse type	12
G	
General safety instructions	8
glas vial	51
Grinding aid.....	40
Grinding chamber inspection	34
Grinding jar	
filling level.....	35
identification	38
inserting.....	27
with safety closure device	41
Grinding jar cleaning.....	38
Grinding jar volume.....	35
Grinding time.....	46
H	
Handling grinding jars	37
Height.....	22
Hoist	
attaching.....	17

I	
Increase of sample volume during grinding process.....	35
Inserting the clamping unit.....	32
Inserting the grinding jar.....	31
Installation.....	11
Installation height.....	12, 18
Installation of the device.....	12, 17, 18
Installation site conditions.....	12
Installing additional equipment.....	51
Interval.....	46
IP40.....	21
L	
Latching bracket wear.....	56
Lid closing.....	21
Lid closing mechanism.....	25
Locking sleeve.....	33
Locking the transport rollers.....	17
LpAeq.....	21
M	
Machine type designation.....	19
Mahlbehälter Auswahl.....	34
Mains connection.....	14
Mains frequency.....	12
Maintenance.....	54
Manual mode.....	46
Manual operation.....	45
Manufacturer's address.....	12
Measurement conditions.....	21
Measurement of noise levels.....	21
Menu structure.....	45
Messages errors.....	50
Moderate or mild injury.....	7
Mounting the transport screws.....	16
N	
Noise levels.....	21
Nominal volume.....	34
Notes on the Manual.....	6
Number of grinding stations.....	20
O	
Opening aid for the clamping unit.....	33
Operating elements and displays.....	25
Operating hours.....	49
Operating modes.....	45
Operating software version.....	49
Operating the device.....	23
Operation.....	15, 42
Overview Table of the Operating Elements and the Display.....	25
Overview table of the parts of the device.....	24
P	
Packaging.....	11
Part number.....	12
Pause time.....	46
Power.....	34
Power failure.....	26, 34
Power failure during grinding.....	34
Power version.....	12
Pressure plate.....	56
Probenmaterialien.....	34
Program 01 to 10.....	45
Programme mode.....	47
property damage.....	7
Protective earth.....	14
Protective equipment.....	21
Q	
Quantity ball.....	35
R	
Rated power.....	20
Receptacle volume.....	20
Regulations for the place of installation.....	14, 20
Releasing the grinding jar clamping mechanism.....	33
Remaining grinding time.....	34
Removing the O-ring.....	38
Repair.....	9
Required floor space.....	22
Residual current protection device.....	14
Return for service and maintenance.....	57
Return goods dispatch note.....	57
Rubber washer.....	56
S	
safety closure device.....	41
Safety locking device	39
Safety notice.....	49
Safety warnings.....	7
Sample quantity.....	35
Sample volumes.....	35
Save parameters.....	47
Selection bar.....	43
Serial number.....	12
serious injury.....	7
Service.....	49
Service Address.....	9
Set language.....	48
shielded cable.....	14
Sintered corundum.....	34
Software update.....	49
Software version of controller.....	49
Software version of display.....	49
Special steel.....	34
Speed.....	46
Speed limits for the adapter.....	53
Stacking the grinding jars.....	36
Stainless steel.....	34
Start in.....	47
Suppressor capacitor wiring.....	14
Symbols in the Display Unit.....	42

T

Target group	19
Target group	8
Technical data	19
Temperature fluctuation and condensed water ..	11
Time	49
Transport	11, 15
Transport lock	
removing from the device	16
Transportation lock	15
Tripping current	14
Tungsten carbide	34
Type plate	12, 14, 20
description	12

U

UKCA marking	12
Ultrafine grinding	35
Unscrewing the transport lock from the transport pallet	16
Use of the device for the intended purpose	19

V

View of the back of the device	24
View of the control panel	25
View of the front of the device	23
View of the menu on the display unit	42
Views of the instrument	23

W

Wear	54
Weight	12, 22
Wet grinding	40
with highly flammable materials	40
Width	22
Workplace-related emissions	21
wrong language	44

Y

Year of production	12
--------------------------	----

Z

Zirconium oxide	34
-----------------------	----

PLANETARY BALL MILL

PM 400 | 20.535.xxxx

EU DECLARATION OF CONFORMITY

We, represented by the undersigned, hereby declare that the above device complies with the following directives and harmonised standards:

Machinery Directive 2006/42/EC

Applied standards, in particular:

DIN EN ISO 12100	Machine Safety - General Design Principles
DIN EN 61010-1	Safety Regulations for Electrical Measurement, Control, Regulation and Laboratory Devices

Electromagnetic compatibility 2014/30/EU (tested at 230 V, 50 Hz)

Applied standards, in particular:

EN 55011	Industrial, scientific and medical equipment - Radio-frequency disturbance characteristics - Limits and methods of measurement
DIN EN 61326-1	Electrical equipment for measurement, control and laboratory use - EMC requirements

Restriction of hazardous substances (RoHS) 2011/65/EU

Authorised person for compilation of the technical documentation:

Julia Kürten (Technical Documentation)

Furthermore, we declare that the relevant technical documentation for the above device has been prepared in accordance with Annex VII Part A of the Machinery Directive and we undertake to submit the documentation to the market surveillance authorities on request.

In the event of a modification of the device not agreed on by Retsch GmbH, as well as the use of non-approved spare parts or accessories, this declaration loses its validity.

Retsch GmbH

Haan, 09/2023



Dr. Frank Janetta, Head of Development





Copyright

© Copyright by
Retsch GmbH
Retsch-Allee 1-5
42781 Haan
Germany